

HD35AP-S / HD35AP-CFR21 release notes

Revision Number	Date	Notes	"IQ/OQ/PQ Guidelines" referenced sections
0.0.4.25	26/02/2021	Minor improvements.	
0.0.4.24	12/02/2021	Added support for HD35AP4GMT.	
0.0.4.22	03/02/2021	Extended the maximum number of measurements managed for each data logger from 12 to 28 (for data loggers that support them, e.g. HD33 series).	
0.0.4.20	11/01/2021	In the „Network options“ section added the possibility to configure the communication with the Delta OHM Cloud and the HDServer1 software.	
0.0.4.18	15/12/2020	Improved TCP server connection type. Added configuration window for LR35... models.	
0.0.4.14	25/11/2020	Added management of new LR35... models. Minor improvements.	
0.0.4.13	11/11/2020	The measurement channels are now displayed with the user-defined labels when viewing data in the database and setting alarms.	
0.0.4.12	05/10/2020	Added possibility to set 12h and 24h as logging interval.	
0.0.4.11	05/05/2020	Added management of “digital input” and “Total energy (kWh)” measurements for data loggers with Modbus RTU input.	
0.0.4.10	24/04/2020	Added HD208 units of measurement management for the japanese market. Added management of new LR35... models.	

Revision Number	Date	Notes	"IQ/OQ/PQ Guidelines" referenced sections
0.0.4.9	06/11/2019	Added the description of the device in alarm in the audit trail. (only HD35AP-CFR21) Added flow units of measurement. Added measurement types for particles counter. Added Modbus RTU Slave management for HD33MT series.	3.7.2 3.11.8
0.0.4.6	22/10/2019	Revision of the user password updating policy in CFR21 environment. (only HD35AP-CFR21) Added new intervals (1, 5 and 10 min) in the periodical transmission to remote servers (FTP, email).	3.7.1 3.7.4 3.11.7 3.11.11
0.0.4.5	14/10/2019	Added new measurement type to support the new HD50CR data logger.	
0.0.4.4	31/07/2019	Added units of measurement management for the japanese market. Added support for centesimal resolution temperature measurements in HD35ED...7P/xTC models. Added Albedo measurement. Speeding up of the alarm window opening. (only HD35AP-CFR21)	3.8.15 3.8.16
0.0.4.2	28/05/2019	Minor changes.	
0.0.4.1	10/04/2019	Added new features for Audit Trail (filtering system) and alarms management. (only HD35AP-CFR21)	3.7.2 3.8.16 3.11.8 3.11.29 3.11.30
0.0.4.0	04/04/2019	HD35AP-S basic software and CFR21 advanced software are now two separate softwares (previously the CFR21 option could be enabled / disabled from the menu inside the single HD35AP-S software).	

Revision Number	Date	Notes	"IQ/OQ/PQ Guidelines" referenced sections
0.0.3.90	26/11/2018	Added daily statistics in the report printing. RH calibration correction for HD208. Minor changes.	
0.0.3.83	12/10/2018	Improvements in the management of the strings of the languages.	