

HD3405.2



HD3405.2 **BENCH-TOP pH METER**

The **HD3405.2** is a bench top instrument for electrochemical measures: **pH**, and temperature. The displayed data can be stored (datalogger) and can be transferred to PC or serial printer thanks to the multi-standard serial port RS232C and USB2.0 and software DeltaLog9 (Vers.2.0 and subsequent ones). The storing and printing parameters can be set from menu.

The HD3405.2 measures pH, redox potential (ORP) in mV. It measures temperature with Pt100 or Pt1000 immersion, penetration or contact probes. The pH electrode calibration can be carried out on one, two or three points and the calibration sequence can be chosen from a list of 13 buffers. The display shows continually the temperature in °C or °F and one of the parameters according to the connected probe type. Printing and storage always include the temperature in °C or °F and one selectable parameter for each probe type.

Other functions include: Max, Min and Avg function, the Auto-HOLD function, the automatic turning off which can also be excluded.

The instruments have IP66 protection degree.

Technical characteristics HD3405.2 pH - mV - °C/°F measurement

Instrument

Dimensions (Length x Width x Height)

Weight

Materials Display

220x120x55mm

460g (complete with batteries)

ABS, rubber

2x41/2 characters plus symbols

visible area: 52x42mm

Operating conditions

Working temperature -5 ... 50°C Storage temperature -25 ... 65°C

Working relative humidity 0 ... 90% RH without condensation

Protection degree

Power

Batteries 3 batteries 1.5V type AA

100 hours with 1800mAh alkaline batteries Autonomy (only batteries) Mains (cod. SWD10) Output mains adapter 100-240Vac/ 12Vdc-1A

Security of memorized data Unlimited

Storage of measured values

Type Quantity

2000 pages of 17 samples each

34,000 sets of measures made up of [pH or mV] and

[°C or °F].

Selectable storage interval 1s, 5s, 10s, 15s, 30s, 1min, 2min, 5min, 10min,

15min, 20min, 30min and 1hour

Time Date and hour Accuracy

Schedule in real time 1min/month max departure

Serial interface RS232C

RS232C electrically isolated Type Baud rate Can be set from 1200 to 38400 baud

Data bit 8 Parity None Stop bit Flow Control Xon/Xoff Serial cable length

Selectable print interval immediate or 1s, 5s, 10s, 15s, 30s, 1min, 2min, 5min,

10min, 15min, 20min, 30min and 1hour

USB Interface

1.1 - 2.0 electrically isolated Type

Connections

Serial interface and USB 8-pole MiniDin connector

Mains adapter (cod. SWD10) 2-pole connector (positive at centre) 12Vdc/1A

Measurement connections Temperature probe input

with SICRAM module or TP47 module

pH/mV input

8-pole male DIN45326 connector

8-pole Mini DIN

female BNC





pH Measurement

Measurement range -2.000...+19.999pH

Resolution 0.01 or 0.001pH selectable from menu

 $\begin{array}{ll} \mbox{Accuracy} & \pm 0.001 \mbox{pH} \pm 1 \mbox{digit} \\ \mbox{Input impedance} & > 10^{12} \mbox{\Omega} \\ \mbox{Calibration error } @25^{\circ} \mbox{C} & \mbox{IOffsetl} > 20 \mbox{mV} \end{array}$

Slope < 50 mV/pH or Slope > 63 mV/pH Sensitivity < 85% or Sensitivity > 106.5%

Automatic / manual -50...+150°C

temperature compensation

mV Measurement

Measurement range -1999.9...+1999.9mV

 Resolution
 0.1mV

 Accuracy
 ±0.1mV ±1digit

 Drift after 1 year
 0.5mV/year

Automatically detected pH standard solutions (@25°C)

1.679pH - 2.000pH - 4.000pH - 4.008pH 4.010pH - 6.860pH - 6.865pH - 7.000pH 7.413pH - 7.648pH - 9.180pH - 9.210pH

10.010pH

Temperature Measurement

 Pt100 measurement range
 -200...+650°C

 Pt1000 measurement range
 -200...+650°C

 Resolution
 0.1°C

 Accuracy
 ±0.1°C ±1digit

 Drift after 1 year
 0.1°C/year

ORDERING CODES

HD3405.2: The kit is composed of: instrument HD3405.2 datalogger, for measurement of pH - redox - temperature, 3 1.5V alkaline batteries, operating manual and DeltaLog9 version 2.0.

pH/mV electrodes, temperature probes, standard reference solutions connection cables for pH electrodes with S7 connector, cables for data download to PC or printer have to be ordered separately.

Accessories

HD2110CSNM: 8-pole connection cable Mini Din - Sub D 9-pole female for RS232C, for connection to PC without USB input.

HD2101/USB: Connection cable USB 2.0 connector type A - 8-pole Mini Din for connection to PC with USB input.

SWD10: Stabilized power supply at 100-240Vac/12 Vdc-1A mains voltage.

HD40.1: Portable, serial input, 24 column thermal printer, 57mm paper width.

HD22.2: Laboratory electrode holder composed of basis plate with incorporated magnetic stirrer, staff and replaceable electrode holder. Height max. 380mm. Powered by benchtop meters of the series HD22... with cable HD22.2.1 (optional) or supplier SWD10 (optional).

HD22.3: Laboratory electrode holder with metal basis plate. Flexible electrode holder for free positioning. For Ø 12mm probes.

TP47: Module for the connection of Pt100 4-wire and Pt1000 2-wire probes to instrument series HD34..., without amplifying electronics.

pH Electrodes

KP 20: Gel pH combined electrode for general use, with S7 screw connector, EPOXY body.

KP 30: Gel pH combined electrode for general use, 1m cable with BNC, EPOXY body .

KP 50: Gel pH combined electrode, porous Teflon ring junction, suitable for emulsions, demineralised water and waste water with S7 screw connector, glass body.

KP 61: 3 diaphragm liquid filled pH combined electrode for wine, milk, cream, etc., S7 screw connector, liquid reference filling, glass body.

KP 62: 1 diaphragm gel pH combined electrode for general use, pure water, varnishes, gel filled, S7 screw connector, glass body.

KP 63: 1 liquid filled pH combined electrode for general use, varnishes, 1m cable with BNC, glass body.

KP 64: Liquid filled pH combined electrode, Teflon ring diaphragm, for wine, varnishes, emulsions, S7 screw connector, glass body.

KP 70: Pointed gel combined pH microelectrode diam. 6 x L=70 mm., with S7 screw connector, EPOXY body, glass tip, open junction for meat and cheese.

KP 80: Pointed gel pH combined electrode, with S7 screw connector, glass body, for cream, milk, viscous material, open junction.

KP100: Flat membrane gel combined pH electrode with S7 screw connector, glass body, for skin, leather, paper.

CP: 1.5m extension cable with BNC/S7 connector for electrode without cable, thread S7.

CP 5: 5m extension cable with BNC/S7 connector for electrode without cable, thread S7.

CP 10: 10m extension cable with BNC/S7 connector for electrode without cable, thread S7.

CP 15: 15m extension cable with BNC/S7 connector for electrode without cable, thread S7.

CE: S7 screw connector for pH electrode.

BNC: female BNC for extension cable

ORP Electrodes

KP90: Redox Platinum electrode, with screw connector S7, electrolyte KCl 3M, body in glass.

KP91: Redox Platinum electrode with 1m cable, GEL fi lled, body in glass.

pH buffer solutions

HD8642: Buffer solution 4.01pH - 200cc. HD8672: Buffer solution 6.86pH - 200cc. HD8692: Buffer solution 9.18pH - 200cc.

Redox buffer solutions

HDR220: Redox buffer solution 220mV 0,5 I. **HDR468:** Redox buffer solution 468mV 0,5 I.

Electrolyte solutions

KCL 3M: 100cc ready for use solution for refilling of the electrodes.

Cleaning and maintenance

HD62PT: Diaphragm cleaning (tiourea in HCl) - 500ml. HD62PP: Protein cleaning (pepsin in HCl) - 500ml. HD62RF: Regeneration (fluorhydric acid) - 100ml. HD62SC: Solution for electrode preservation - 500ml.

Temperature probes equipped with SICRAM module

TP472I: Wire wound Pt100 sensor, immersion probe. Stem \emptyset 3 mm, length 300 mm. Cable length 2 m.

TP472I.0: Thin film Pt100 sensor, immersion probe. Stem Ø 3 mm, length 230 mm. Cable length 2 m.

TP473P.I: Wire wound Pt100 sensor, penetration probe. Stem Ø 4mm, length 150 mm. Cable length 2 m.

TP473P.O: Thin film Pt100 sensor, penetration probe. Stem Ø 4mm, length 150 mm. Cable length 2 m.

TP474C.I: Wire wound Pt100 sensor, contact probe. Stem Ø 4mm, length 230mm, contact surface Ø 5mm. Cable length 2 m.

TP474C.0: Thin film Pt100 sensor, contact probe. Stem Ø 4mm, length 230mm, contact surface Ø 5mm. Cable length 2 m.

TP475A.0:, Thin film Pt100 sensor, air probe. Stem Ø 4mm, length 230mm. Cable length 2 m. TP472I.5: Thin film Pt100 sensor, penetration probe. Stem Ø 6mm, length 500 mm. Cable length 2 m.

TP472I.10: Thin film Pt100 sensor, penetration probe. Stem Ø 6mm, length 1000mm. Cable length 2 m.

TP49A.0: Thin film Pt100 sensor, immersion probe. Stem Ø 2,7mm, length 150mm. Cable length 2 m. Aluminium handle

TP49AC.0: Thin film Pt100 sensor, contact probe. Stem Ø 4mm, length 150mm. Cable length 2 m. Aluminium handle

TP49AP.0: Thin film Pt100 sensor, penetration probe. Stem Ø 2,7mm, length 150mm. Cable length 2 m. Aluminium handle

TP875.I: Wire wound Pt100 sensor, 150mm diameter globe-thermometer equipped with handle. Cable length 2 m.

TP876.I: Wire wound Pt100 sensor, 50mm diameter globe-thermometer equipped with handle. Cable length 2 m.

TP87.0: Thin film Pt100 sensor, immersion probe. Stem Ø 3 mm, length 70 mm. Cable length 2 m.

TP878.0: Thin film Pt100 sensor, contact probe for solar panels. Cable length 2 m.

TP878.1.0: Thin film Pt100 sensor, contact probe for solar panels. Cable length 5 m.

TP879.0: Thin film Pt100 sensor, penetration probe for compost. Stem Ø 8 mm, length 1000 mm. Cable length 2 m.

Temperature probes without SICRAM module

TP47.100.0: Thin film Pt100 sensor, immersion probe. Stem Ø 3mm, length 230mm. Connection cable 4 wires with connector, length 2 m.

TP47.1000.0: Thin film Pt1000 sensor, immersion probe. Probe's Stem Ø 3mm, length 230mm. Connection cable 4 wires with connector, length 2 m.

TP47: Connector for Pt100 4-wire and Pt1000 2-wire probes without SICRAM module.

TP87.100.0: Thin film Pt100 sensor, immersion probe. Stem Ø 3mm, length 70mm. 4-wires connection cable with connector, length 1 m.

TP87.1000.0: Thin film Pt1000 sensor, immersion probe. Stem Ø 3mm, length 70mm. 2-wires connection cable with connector, length 1 m.



