

# English

# **Operating manual**

Conductivity meters / pH meters / thermometers **HD2156.1 – HD2156.2** 



Members of GHM GROUP:

GREISINGER HONSBERG *Martens Selta*cers VAL.CO

www.deltaohm.com Keep for future reference.

INTRODUCTION	
KEYBOARD AND MENU DESCRIPTION	8
PH MEASUREMENT	16
Automatic or manual pH compensation PH ELECTRODE CALIBRATION Temperature characteristics of Delta OHM pH standard solutions	17
CONDUCTIVITY MEASUREMENT	21
Automatic or manual conductivity compensation Calibration of conductivity Table of standard solution at 147μS/cm, 1413μS/cm, 12.88mS/cm and 111800mS/cm	23
DIRECT INPUT INTO Pt100 AND Pt1000 TEMPERATURE PROBES How to measure Instructions to connect the TP47 module for conductivity/temperature combined probes, 4-wire Pt100 probes, and 2-wire Pt1000 probes	26
WARNINGS AND OPERATING INSTRUCTIONS	30
Notes on the use of the pH electrodes Notes on conductivity probes INSTRUMENT SIGNALS AND FAULTS	31
LOW BATTERY WARNING AND BATTERY REPLACEMENT	35
INSTRUMENT STORAGE	
SERIAL INTERFACE AND USB	37
STORING AND TRANSFERRING DATA TO A PERSONAL COMPUTER	
The <i>Logging</i> function - only for HD2156.2 Clearing the memory - only for HD2156.2 The <i>PRINT</i> function	
CONNECTION TO A PC	41
CONNECTION TO THE RS232C SERIAL PORT OF THE INSTRUMENT CONNECTION TO THE USB 2.0 PORT OF THE INSTRUMENT - ONLY FOR HD2156.2 NOTES ABOUT WORKING AND OPERATIVE SAFETY	41
INSTRUMENT TECHNICAL CHARACTERISTICS         TECHNICAL DATA OF PROBES         pH ELECTRODES         2 AND 4-ELECTRODE CONDUCTIVITY PROBES         Pt100 4-WIRE PROBES AND Pt1000 2-WIRE PROBES COMPLETE WITH TP47 MODULE	
ORDER CODES	

# CONTENTS

# INTRODUCTION

The **HD2156.1** and **HD2156.2** are portable instruments with a large LCD display. They measure pH, mV, redox potential (ORP), conductivity, liquid resistivity, total dissolved solids (TDS) and salinity using combined 4-ring and 2-ring conductivity/temperature probes. Temperature only is measured by Pt100 or Pt1000 immersion, penetration or contact probes.

The pH electrode calibration, as well as manual, can be carried out automatically on one, two or three points and the calibration sequence can be chosen from a list of 13 buffers. The probe calibration can be performed automatically in one or more of the 147  $\mu$ S/cm, 1413  $\mu$ S/cm, 12880  $\mu$ S/cm or 111800  $\mu$ S/cm conductivity calibration solutions.

The HD2156.2 instrument is a **datalogger**. It memorizes up to 20,000 sets of three measurements composed of pH or mV, conductivity or resistivity or TDS or salinity and temperature: these data can be transferred from the instrument connected to a PC via the RS232C serial port or USB 2.0 port. The storing interval, printing, and baud rate can be configured using the menu.

The HD2156.1 and HD2156.2 models are fitted with an RS232C serial port and can transfer the acquired measurements to a PC or to a portable printer in real time. Printing and, for the model HD2156.2, storing always include three parameters: temperature – pH or mV –  $\chi$  or  $\Omega$  or TDS or g/l.

The *Max*, *Min* and *Avg* function calculates the maximum, minimum or average values. Other functions include: the Auto-HOLD function and the automatic turning off which can also be disabled.

The instruments have IP66 protection degree.


This manual describes the HD2156.1 and HD2156.2 models: if not otherwise specified, the description is intended to be applicable to both models.



Conductivity meter – pH meter – Thermometer HD2156.1

# HD2156.1

- 1. BNC connector for the pH/mV electrode.
- 2. 8-pole DIN45326 connector, input for combined 4-ring or 2-ring conductivity/temperature probes, for direct 4-wire Pt100 temperature probes and 2-wire Pt1000 probes complete with TP47 module.
- 3. External auxiliary power supply connector input (positive at centre).
- 4. Battery symbol: displays the battery charge level.
- 5. Function indicators.
- 6. Secondary display line.
- 7. CAL/▲ key: during normal operation starts the conductivity probe or pH electrode calibration; in the menu, increases the current value.
- 8. **FUNC/ENTER** key: during normal operation displays the maximum (MAX), the minimum (MIN) and the average (AVG) of current measurements; in the menu, confirms the current selection. This updates the display measurement when the *Auto-Hold* function is enabled.
- 9. **pH/mV-ESC** key: changes the main variable measurement between pH and mV. In the menu, confirms the current selection.
- 10. SERIAL key: starts and ends the data transfer to the serial communication port.
- 11. **MENU** key: allows access to and exit from the menu.
- 12. **X-Ω-TDS/ESC** key: changes the main variable measurement between conductivity, resistivity, total dissolved solids (TDS) and salinity. In the menu, cancels the operation in progress without making changes.
- 13. °C/°F key: when the probe is not connected, allows manual modification of the temperature. When the probe is connected, if this key is pressed twice, it changes the unit of temperature measurement from degrees Celsius to Fahrenheit.
- 14. **ON-OFF/AUTO-OFF** key: turns the instrument on and off; when pressed together with the CAL key, disables the automatic turn off.
- 15. pH electrode efficiency indicators.
- 16. MAX, MIN and AVG symbols.
- 17. Main display line.
- 18. Line for symbols and comments.
- 19. 8-pole MiniDin connector for RS232C. For the connection to PC (with cable HD2110CSNM or C206) or printer (with cable HD2110CSNM).

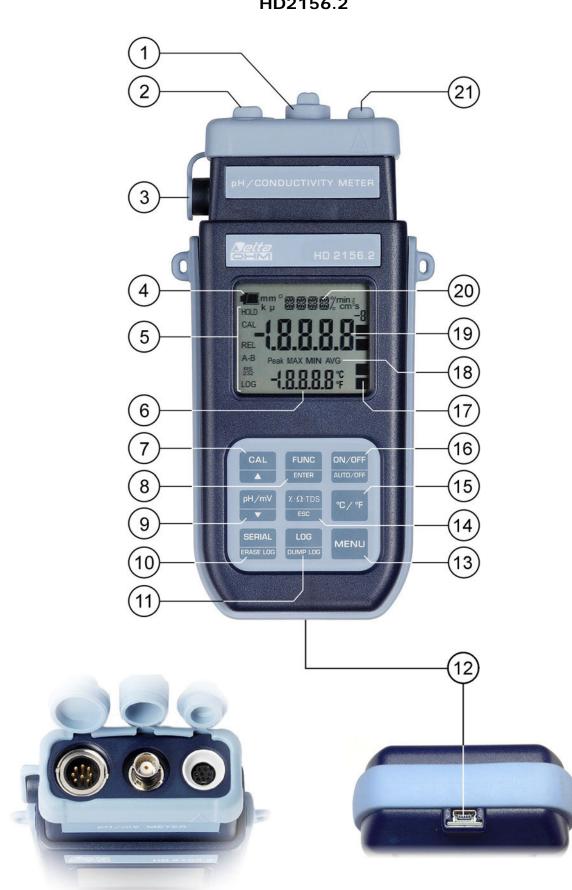

Conductivity meter – pH meter – Thermometer HD2156.2

# HD2156.2

- 1. BNC connector for the pH/mV electrode.
- 2. 8-pole DIN45326 connector, input for combined 4-ring or 2-ring conductivity/temperature probes, for direct 4-wire Pt100 temperature probes and 2-wire Pt1000 probes complete with TP47 module.
- 3. External auxiliary power supply connector input (positive at centre).
- 4. Battery symbol: displays the battery charge level.
- 5. Function indicators.
- 6. Secondary display line.
- 7. CAL/▲ key: during normal operation starts the conductivity probe or pH electrode calibration; in the menu, increases the current value.
- 8. **FUNC/ENTER** key: during normal operation displays the maximum (MAX), the minimum (MIN) and the average (AVG) of current measurements; in the menu, confirms the current selection. This updates the display measurement when the *Auto-Hold* function is enabled.
- 9. **pH/mV-ESC** key: changes the main variable measurement between pH and mV. In the menu, confirms the current selection.
- 10. **SERIAL/ERASE LOG** key: starts and ends the data transfer to the serial communication port. In the menu, clears the data contained in the instrument's memory.
- 11. **LOG/DUMP LOG** key: during normal operation, starts and ends the saving of the data in the internal memory; in the menu, starts the data transfer from the instrument's memory to the PC.
- 12. Mini-USB type B connector for USB 2.0. For the connection to PC (with cable CP23).
- 13. MENU key: allows access to and exit from the menu.
- 14. **X-Ω-TDS/ESC** key: changes the main variable measurement between conductivity, resistivity, total dissolved solids (TDS) and salinity. In the menu, cancels the operation in progress without making changes.
- 15. °C/°F key: when the probe is not connected, allows manual modification of the temperature. When the probe is connected, if this key is pressed twice, it changes the unit of temperature measurement from degrees Celsius to Fahrenheit.
- 16. **ON-OFF/AUTO-OFF** key: turns the instrument on and off; when pressed together with the CAL key, disables the automatic turn off.
- 17. pH electrode efficiency indicators.
- 18. MAX, MIN and AVG symbols.
- 19. Main display line.
- 20. Line for symbols and comments.
- 21. 8-pole MiniDin connector for RS232C. For the connection to PC (with cable HD2110CSNM or C206) or printer (with cable HD2110CSNM).

#### **KEYBOARD AND MENU DESCRIPTION**

#### Foreword

The instrument keyboard is composed of single-function keys, like the MENU key, and double-function keys such as the ON-OFF/Auto-OFF key.

In the double-keys, the function in the upper part is the "main function", while the one in the bottom part is the "secondary function". When the instrument is in standard measurement mode, the main function is active. In the menu or in conjunction with the FUNC key, the secondary function is enabled.

The pressing of a key is accompanied by a short confirmation beep: a longer beep sounds if the wrong key is pressed.

Each key specific function is described in detail below.

#### The following can be seen on the display at the same time:


- 1. pH (or mV) and temperature
- 2. conductivity (or resistivity or total dissolved solids or salinity) and temperature.

It is not possible to see pH (or mV) and conductivity at the same time: one measurement or the other must be selected using the "pH/mV -  $\star$ " and "X- $\Omega$ -TDS/ESC" keys. In direct print and when logging three quantities are captured at the same time: pH/mV, conductivity and temperature.



# ON-OFF/Auto-OFF key

The instrument is turned on and off using the ON/OFF key. Turning on enables all display segments for a few seconds, starts an auto-test, and displays the current values of cell constant (CELL) and temperature coefficient  $\alpha$  (ALPH). Finally, the instrument is set in standard working mode, displaying the measurement which remained enabled at the last turning off in the main line, and the temperature in the secondary line.

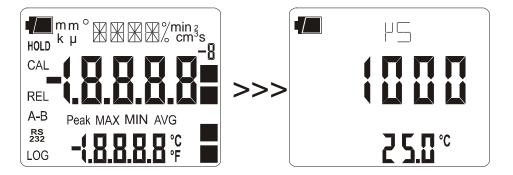

During turning on, should no probes be connected, the last manually-set temperature appears in the secondary line. The unit of measurement symbol (°C or °F) starts blinking, and a letter "m" meaning "manual" appears next to the battery symbol. The probe's data are captured upon turning the instrument on: if the ERR message appears in the secondary line, it is necessary to turn the instrument off and then on again.

Replace the probes when the instrument is off.



The instrument has an *AutoPowerOff* function that automatically turns the instrument off after about 8 minutes if no key is pressed during the intervening time. The *AutoPowerOff* function can be disabled by holding the CAL/ $\checkmark$  key pressed down when turning the instrument on: the battery symbol will blink to remind the user that the instrument can only be turned off by pressing the <ON/OFF> key.

The automatic turning off function is disabled when external power is used. On the other hand, it cannot be disabled when the batteries are discharged.



# FUNC/ENTER key

During normal measurement this enables the display and logging of the maximum (MAX), minimum (MIN) and average (AVG) value of the pH, mV, conductivity, liquid resistivity, total dissolved solids, salinity and temperature measurements, updating them with the acquisition of new samples. The acquisition frequency is once a second.

To switch from pH to mV, press the "pH/mV- $\bullet$ " key. Use the  $\chi$ - $\Omega$ -TDS/ESC key to switch from conductivity to liquid resistivity, to total dissolved solids or to salinity.

In the menu, the ENTER key confirms the current parameter and then goes to the next one.

The MAX, MIN and AVG measurements remain in the memory until the instrument is on, even after exiting the calculation function. To reset the previous values and restart with a new measurement session, press FUNC until the message "FUNC CLR" appears, then use the arrows to select YES and confirm using ENTER.

According to settings in the "RCD Mode" menu item, the maximum, minimum and average indications have different meanings: please see the description of this MENU key below.

Attention: the data captured using the Record function cannot be transferred to the PC.




It increases the current parameter when used in the menu; when used in measurement mode, it starts the displayed variable (pH or conductivity) calibration (please see the chapter dedicated to pH, page 17, and conductivity, page 23, calibration).



When a probe with temperature sensor is connected, the key changes the unit of measurement from degrees Celsius to Fahrenheit. The measured temperature value is used to compensate the pH or conductivity measurement.

If the probe is not present, the compensation temperature must be entered manually: to manually change the value shown in the display lower line, press °C/°F once. The temperature indicated starts blinking. While the display is blinking, it is possible to change the compensation temperature using the arrows ( $\blacktriangle$  and  $\checkmark$ ). Confirm using ENTER. The display stops blinking, and the temperature displayed is used for compensation.

If the temperature probe is not present, to change the unit of measurement between  $^{\circ}C$  and  $^{\circ}F$ , it is necessary to press **twice** the  $^{\circ}C/^{\circ}F$  key.

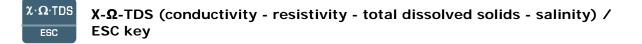

Changes the main variable measurement between conductivity, resistivity, total dissolved solids (TDS) and salinity. The selected parameter is used for display, printing, and logging. In the menu, the key clears or cancels the active function (ESC).

#### Auto-HOLD function

The instrument has an Auto-Hold function, which can be set in the MENU, that "freezes" the measurement automatically when it has been stable (**within 1 digit**) for over 10 seconds: the message HOLD is displayed.

To perform a new measurement, it is necessary to press the FUNC/ENTER key.

The HOLD message starts blinking, while the display follows the actual measurement trend, until it stabilizes again and the HOLD message remains still.



pH/mV - <del>v</del> key

Changes the main variable measurement between pH and mV. The selected parameter is used for display, printing, and logging.

#### Auto-HOLD function

The instrument has an Auto-Hold function, which can be set in the MENU, that "freezes" the measurement automatically when it has been stable (within 1 mV) for over 10 seconds: the message HOLD is displayed.

To perform a new measurement, it is necessary to press the FUNC/ENTER key.

The HOLD message starts blinking, while the display follows the actual measurement trend, until it stabilizes again and the HOLD message remains still.

NOTE: when the Auto-Hold function is enabled, the measurement in mV is disabled. To restore the display in mV, disable the Auto-Hold function in the menu.

When used in the menu, it decreases the current variable value.



The first menu item is accessed by initially pressing on the MENU key; press ENTER to go to the following items. To modify the item displayed, use the arrow keys ( $\checkmark$  and  $\checkmark$ ). The current value is confirmed by pressing the ENTER key and the display moves on to the next parameter. If pressing ESC, the setting is cancelled.

To exit the menu, press the MENU key at any time.

The menu items are listed in this order:


1) Management of memorized data (only HD2156.2): the message "LOG\_DUMP\_or\_ERAS" (Transfer data or erase) is scrolled in the comment line. The center figure reports the number of free memory pages (FREE). All memory data are permanently erased by pressing SERIAL/EraseLOG. By pressing LOG/DumpLOG, the data transfer of the logged data on the serial port is started: the "BAUD-RATE" must have previously been set to the maximum value (please see the menu items described below and the paragraph "STORING AND TRANSFERRING DATA TO A PERSONAL COMPUTER" on page 39).

2) Identifier of the sample being measured: it is an automatically increased progressive number associated with the single PRINT function (print interval set to 0) for the printing of labels. The index appears in the single sample printing together with date, time, conductivity (liquid resistivity, total dissolved solids or salinity) and temperature or pH (or mV) measured values. This menu item allows the value of the first sample to be set: each time the PRINT key is pressed, the identification ID in the printing is increased by 1 allowing progressive measurement of all measured samples. If the Auto-Hold function, described below in this chapter, is enabled, the print time interval is forced to zero. Pressing SERIAL only causes the print to occur when the measurement has stabilized (HOLD symbol still). Later, it is possible to repeat the print at will, but while the HOLD mode is on, the sample identifier number is not increased. This is useful when more labels must be printed with the same identification code without increasing the code each time.

The message "SMPL ID UNT=RSET SER=PRNT" is scrolled in the comment line: using the arrows ( $\blacktriangle$  and  $\checkmark$ ) the currently measured sample identifier value can be changed. By holding the °C/°F (Unit) key down the proposed number is rapidly set to zero. The instrument's heading information will be printed using the SERIAL key.

3) **AUTO-HOLD function**: the instrument normally operates in **continuous view** mode (default setting). In this mode the displayed measurement is updated every second. If the Auto-Hold function is enabled, the instrument performs the measurement and when it stabilizes it goes in HOLD mode. To update the display indication, press FUNC/ENTER. The Auto-Hold function is applied to the pH measurement, conductivity and all resulting measurements.

In the following figure you can see an example of the measurement process with the Auto-Hold function enabled. A probe is immersed into a liquid at conductivity  $\chi_1$  and, to perform the measurement, the FUNC/ENTER key is pressed: The conductivity measurement raises progressively reaching the final value. The HOLD symbol blinks. In the stretch indicated by 1, the measurement remains stable for 10 seconds, within two digits: at the end of this interval (point 2), the instrument goes into HOLD mode, presenting the final stable value.

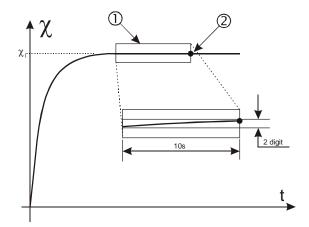

4) Print and log interval: sets the interval in seconds between two loggings or data transfers to the serial port. The interval can be set at 0, 1s, 5s, 10s, 15s, 30s, 60s (1min), 120s (2min), 300s (5min), 600s (10min), 900s (15min), 1200s (20min), 1800s (30min) and 3600s (1hour). If the value 0 is set, SERIAL works on command: the sending of data to the serial port is performed each time the key is pressed. Recording (LOG) is performed with one second intervals even if the interval is set to 0. With an interval from 1 to 3600s, continuous data transfer is started when the SERIAL key is pressed. To end the recording (LOG) and continuous data transfer operations (SERIAL with an interval greater than 0), press the same key again.

5) Sleep\_Mode\_LOG (Automatic turning off during recording) (only HD2156.2): this function controls the instrument's automatic turning off during logging, occurring between the capture of a sample and the next one. When the interval is lower than 60 seconds, the instrument will always remain on. With intervals greater than or equal to 60 seconds, it is possible to turn off the instrument between loggings: it will turn on at the moment of sampling and will turn off immediately afterwards, thus increasing the battery life. Using the arrows select YES and confirm using ENTER in order to enable the automatic turning off, select NO and confirm to disable it and keep the instrument on continuously.

Note: even if **Sleep\_Mode\_LOG=YES** is selected, the instrument does not turn off for less than one minute intervals.

- 6) **LAST CAL pH m/d h/m (Last pH calibration)**: the display shows the month and day (m/d) in the main line, and the hour and minutes (h/m) in the secondary line of the last pH electrode calibration. This menu item cannot be modified. The calibration year is not displayed.
- 7) **BUFR\_1\_pH (First pH buffer)**: selects the value of the first buffer for the pH electrode calibration. The preset buffers are compensated for temperature. The ATC indication is displayed in the lower line. The USER buffer, defined by the user, is not compensated for temperature: to change its value, go to the "SET USER BUFR pH" step. Please see the paragraph dedicated to pH calibration on page 17.
- 8) **BUFR\_2\_pH (Second pH buffer)**: selects the value of the second buffer for the pH electrode calibration. No buffer solution values too close (less than 2pH) to the first buffer are proposed. This is also true for the USER buffer, defined by the user at the "SET USER BUFR pH" step: if, for example BUFR\_1=6.860 and USER=5.000, the BUFR\_2 USER does not appear among the values because it is too close to BUFR\_1. The selection NIL (=no buffer) disables the current buffer. The USER buffer, defined by the user, is not compensated for temperature. Please see the paragraph dedicated to pH calibration on page 17.
- 9) BUFR\_3\_pH (Third pH buffer): selects the value of the third buffer for the pH electrode calibration. No buffer solution values too close (less than 2pH) to the first two buffers are proposed. This is also true for the USER buffer, defined by the user at the "SET USER BUFR pH" step: The selection NIL (=no buffer) disables the current buffer. The USER buffer is not compensated for temperature. Please see the paragraph dedicated to pH calibration on page 17.
- 10) **SET USER BUFR pH (Set user pH buffer)**: this is a buffer whose value can be defined by the user using the arrows; all values from 0 to 14pH are available. This buffer is not compensated for temperature, so the buffer value must be set at the actual solution temperature. As an alternative, the correct value according to temperature can be set in the calibration phase.
- 11) **K CELL (Conductivity cell constant):** sets the conductivity probe's cell constant nominal value. The values 0.01, 0.1, 0.7, 1.0 and 10 cm<sup>-1</sup> (with tolerances from –30% to +50% of nominal value) are admitted. The cell constant must be inserted before starting the probe calibration. An ERR signal is generated if the cell constant's actual value exceeds the limits –30% or +50% of nominal value. In this case it is necessary to check that the value set is correct, that the calibration solutions are in good state, and then proceed with a new calibration.

The cell constant change entails resetting the calibration date: a new calibration updates the calibration date.

12) **LAST CAL K COND m/d h/m (Last conductivity calibration)**: the display shows the month and day (m/d) in the main line, and the hour and minutes (h/m) in the secondary line of the previous calibration of the conductivity probe. This menu item cannot be modified. The calibration year is not displayed. The cell constant change using the K CELL\_COND parameter, resets the date.

- 13) ALPH\_T\_COND (Temperature coefficient α<sub>T</sub>): the temperature coefficient α<sub>T</sub> is the percentage measurement of the conductivity variation according to temperature and is expressed in %/°C (or %/°F). The admitted values vary from 0.00 to 4.00%/°C. Use the arrows (▲ and ◄) to set the desired coefficient α<sub>T</sub>, and confirm with ENTER.
- 14) **REF\_TEMP\_COND (Conductivity reference temperature)**: it indicates the temperature to which the displayed conductivity value is standardized and can be equal to 20°C or 25°C. Using the arrows (▲ and ◄), set the desired value and confirm with ENTER.
- 15) **TDS COND (Conversion factor X/TDS)**: it represents the ratio between the measured conductivity value and the total quantity of dissolved solids in the solution, expressed in mg/l (ppm) or g/l (ppt). This conversion factor depends on the nature of the salts present in the solution: in the field of water quality treatment and control, where the main component is CaCO<sub>3</sub> (Calcium Carbonate), a value of 0.5 is usually used. For agriculture water, for fertilizers preparation, and in hydroponics, a factor of about 0.7 is used. Using the arrows (▲ and ▼), set the desired value, selecting it in the 0.4...0.8 range, and confirm with ENTER.
- 16) **RCD MODE (Record mode)**: the instrument captures a pH, a conductivity and a temperature value every second. If the RCD MODE parameter is set to "**conductivity**", the maximum (MAX) and minimum (MIN) values displayed using FUNC/ENTER refer to conductivity: the indicated temperature and pH are those measured at the maximum and minimum conductivity and are not the maximum and minimum temperature and pH. Similarly, if the RCD MODE parameter is set to "**pH**", the maximum and minimum values displayed using FUNC/ENTER refer to pH: the indicated temperature and conductivity are those measured at the maximum and minimum pH and are not the maximum and minimum temperature and conductivity.

If the RCD MODE parameter is set to "**tp**" (=temperature), the maximum and minimum values displayed using FUNC/ENTER refer to temperature: the indicated conductivity and pH are those measured at the maximum and minimum temperature and are not the maximum and minimum conductivity and/or pH.

Finally, if the RCD MODE parameter is set to "**Indep**" (=independent), the maximum and minimum values displayed using FUNC/ENTER are independent: the indicated pH, conductivity and temperature are the maximum and minimum measured values but are not necessarily referred to the same measurement moment (factory default).

- 17) **Probe type**: the message "PRBE\_TYPE" is scrolled in the comment line. The main line in the center of the display shows the type of temperature probe connected to the instrument. Conductivity/temperature combined probes with Pt100 or Pt1000 sensor, or temperature only probes can be connected to the input.
  - 4-wire Pt100 using the TP47 module
  - 2 or 4-wire Pt1000 using the TP47 module

**Upon being turned on,** the instrument automatically detects the temperature probes: the *Probe Type* menu item is configured by the instrument and cannot be modified by the user. If no temperature probe or combined probe with temperature sensor is connected, the instrument selects the Pt1000 sensor.

- 18) **YEAR**: to set the current year. Use the arrows to modify this parameter and confirm using ENTER.
- 19) **MNTH (month)**: to set the current month. Use the arrows to modify this parameter and confirm using ENTER.
- 20) **DAY**: to set the current day. Use the arrows to modify this parameter and confirm using ENTER.
- 21) **HOUR**: to set the current hour. Use the arrows to modify this parameter and confirm using ENTER.

- 22) **MIN**: to set the current minutes. In order to correctly synchronize the minute, it is possible to reset the seconds by pressing the °C/°F key. Use the arrows to set the current minute plus one, and as soon as that minute is reached press °C/°F: this synchronizes the time to the second. Press ENTER to go onto the next item.
- 23) BAUD\_RATE: indicates the frequency used for the serial communication with the PC. Values from 1200 to 38400 baud. Use the arrows to modify this parameter and confirm using ENTER. The communication between instrument and PC (or serial port printer) only works if the instrument and PC baud rates are the same. If the USB connection is used this parameter value is automatically set (please see the details on page 39).



# LOG/DumpLOG key - only HD2156.2

In measurement mode, this function starts and stops the logging of a data block to be saved in the instrument's internal memory. The data logging frequency is set in the "**Print and log in-terval**" menu parameter. The data logged between a start and subsequent stop represent a block.

When the logging function is on, the LOG indication is displayed, the battery symbol blinks and a beep is issued each time a logging occurs; **the battery symbol does not appear when using an external power supply**.

To end the logging, press LOG.


# If the Auto-HOLD function is enabled (please see the menu), the data logging is disabled.

The HD2156.2 can turn off during logging between one capture and the next: the function is controlled by the **Sleep\_Mode\_LOG** parameter. When the logging interval is less than one minute, the logging instrument remains on; with an interval of at least one minute, it turns off between one capture and the next if the parameter **Sleep\_Mode\_LOG**=YES.



When the LOG key is pressed after the MENU key, the transfer of the logged data on the serial port is started.

Please see the paragraph dedicated to data transfer on page 39.

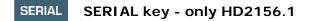



SERIAL/EraseLOG key - only HD2156.2

In measurement mode, this function starts and stops the data transfer to the RS232C serial output.

According to the settings entered in the **Print and log interval** menu item, a single sample can be printed if **Print and log interval**=0 or a continuous indefinite printing of the measured data can be set up if **Print and log interval**=1...3600.

The printing operation is accompanied by the display of the RS232 symbol and the blinking of the battery symbol; when using an external power supply the battery symbol does not appear.

Press SERIAL to end the continuous printing.

Before starting the printing with SERIAL, set the baud rate. To do so, select the **Baud Rate** menu item and select the maximum value equal to 38400 baud by using the arrows. Confirm by pressing ENTER.

The DeltaLog9 software for PC will automatically set the baud rate value during connection. If you are using a different program than DeltaLog9, be sure the baud rate is the same for both the instrument and the PC: the communication will only work in this way.



When pressed after the MENU key, the SERIAL key **permanently** erases all the data contained in the instrument's memory.

#### **pH** MEASUREMENT

The instruments work with pH measurement electrodes, redox potential measurement electrodes (ORP), and specific ion electrodes. The pH measurement must always be accompanied by a temperature measurement. Probes with Pt100 or Pt1000 sensor may be used for measuring temperature or for the automatic compensation of the Nernst coefficient with the pH electrode.

The pH or mV indication is displayed in the main line by pressing the "pH/mV" key; the secondary line shows the temperature. If no temperature probe with temperature sensor is connected, the secondary line shows the manual temperature.

#### The electrode for pH measurement

The electrode for pH measurement, generally in glass, generates an electrical signal proportional to the pH according to Nernst law. Of this signal the following aspects are considered:

*Zero point*: The pH where the electrode generates a potential of 0 mV. In most electrodes, this value is found at about 7pH.

*Offset or Asymmetry Potential*: mV generated by an electrode when immersed in a standard solution at 7pH. Generally oscillates between ± 20 mV.

*Slope*: response of the electrode expressed in mV per pH units. The theoretical electrode slope at 25°C is 59.16 mV/pH. When the electrode is new the slope is close to the theoretical value.

*Sensitivity*: it is the electrode's slope expression in relative terms. It is obtained by dividing the actual value of the slope by the theoretical value, and is expressed as a %. The asymmetry potential and the slope vary in time with the use of the electrode, which necessitates regular calibration.

The pH electrodes must be calibrated using the standard solutions (see the pH calibration chapter below). The ORP and specific ion electrodes do not need calibration as their absolute voltage is measured. The standard redox solutions are only used to check the quality of a redox electrode.


There is no user temperature calibration.

#### The probes are detected during turn on, and this cannot be performed when the instrument is already on, therefore if a probe is connected and the instrument is on, it is necessary to turn it off and on.

#### AUTOMATIC OR MANUAL pH COMPENSATION

In a correct measurement of pH, the results need to be expressed together with the temperature value at which the reading is performed.

The electrode slope varies according to the temperature in a known mode according to Nernst law: e.g., a 1 pH variation, that at 25 °C means 59.16 mV, at 100 °C means 74.04 mV.

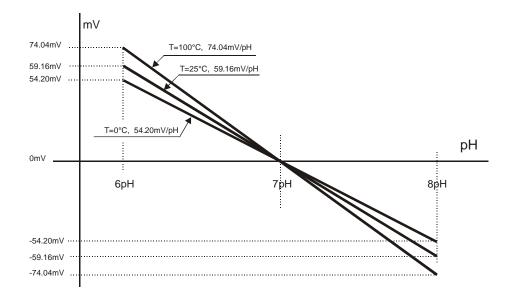

When a temperature probe is present, the instrument automatically applies the ATC automatic temperature compensation function. To disable it, the temperature probe must be disconnected.

If the temperature probe is not present, and the correct value is not entered manually, the extent of the error committed in pH measurement is proportional to temperature and pH value itself.

In absence of the temperature probe, the lower display shows the manually set compensation temperature (default=25 °C).

To point this condition out, the °C or °F symbol blinks intermittently near the temperature value. Moreover, on the main display an "m" (manual) is turned on near the battery symbol (if present). On the print-outs, the **MT** indication is printed. Instead, if the probe is present, the **AT** symbol is printed.



To manually change the compensation temperature press °C/°F once: the indicated temperature value starts blinking. Select the desired temperature value by using the arrows and confirm with ENTER. The display stops blinking, and the temperature displayed is used for compensation.

To change the unit of measurement between °C and °F, it is necessary to press twice the °C/°F key.

# PH ELECTRODE CALIBRATION

The electrode calibration is used to compensate the zero potential and slope deviations to which the electrode is subject with time.

The calibration frequency depends on the accuracy desired by the user and by the effects that the measured sample have on the electrode. Generally, we recommend daily calibration, but it is the user's responsibility from personal experience, to establish the most appropriate frequency.

The calibration may be carried out using 1, 2 or 3 points. When using 1 point, the electrode offset is corrected, with 2 points the offset and the gain is corrected; finally, in case a three points calibration is carried out two offsets and two slopes are calculated, **the second point is the one in which the zero is performed**.

The instrument has a memory of 13 buffers with relevant temperature compensation tables (ATC) plus an "User" buffer, not compensated. The three buffers can be selected by using the **BUFR\_1\_pH, BUFR\_2\_pH, BUFR\_3\_pH** menu items. Usually one for the acid, one for the neutral, and one for the alkaline band will be selected:

@25 C					
BUFR_1_pH (NEUTRAL)	6.860	6.865	7.000	7.413	7.648
BUFR_2_pH (ACID)	1.679	2.000	4.000	4.008	4.010
BUFR_3_pH (ALKALINE)	9.180	9.210	10.010		

The buffers in bold are predefined in the factory, these are the "DeltaOHM" ones.

If electrode calibration was not carried out on the instrument, or the batteries have been changed or the last calibration failed, the display blinks the **CAL** message.

## Calibration procedure

താടംറ

- 1. Display the measurement in pH by pressing the "pH/mV" key.
- 2. Insert the temperature probe and the electrode in the solution of a calibration buffer. If no temperature probe is available, use another thermometer and enter the value manually as indicated in the paragraph "Automatic or manual pH compensation".
- 3. The electrode calibration mode is started by pressing CAL.
- 4. Among the three prepared buffers, the instrument automatically detects the closest to the pH value being read and flags it up on the lower display. If the buffer is not detected the first buffer BUFR\_1\_pH is proposed.
- 5. At this point the CAL symbol is not blinking on the display and the following is shown from the upper side downward:
  - a) the chosen buffer nominal value at 25 °C (scrolling indication)
  - b) the pH measurement value with the current calibration
  - c) the temperature compensated buffer value
- 6. The detected and temperature compensated buffer value, shown in the lower line, can be modified using the arrows.
- 7. To proceed with the calibration press ENTER. The electrode offset mV value (OFFS) is shown on the display for few seconds.
- 8. After a few seconds the instrument shows the measurement again, corrected according to the new calibration, **but remains in calibration mode**. Pressed repeatedly the ENTER key allows the calibration on the point to be repeated, for example, in order to obtain a more stable value.
- 9. To end the electrode calibration, press "ESC" to exit calibration, or continue the calibration for the second point.
- 10. Extract the electrode from the buffer, clean it carefully, and insert it in the following buffer.
- 11. Press the MENU key.
- 12. The instrument displays the value detected on the new buffer: continue by repeating the steps from point 4.

# NOTES:

- After calibration, the instrument displays an electrode quality indication:
  - No signal: electrode functioning.
  - 1 small square blinking on the lower right: electrode almost exhausted.
  - 2 small squares blinking on the lower right: electrode exhausted to be replaced.
- The 3 point calibration must always be carried out according to the fixed sequence: NEUTRAL>>ACID>>BASIC. The basic buffer must be the last in the sequence.
- Without having pressed ENTER at all, the calibration is interrupted by pressing ESC; the previous values will continue to be used.
- The buffers are always presented in the sequence set in the menu with the BUFR\_1-BUFR\_2-BUFR\_3 parameters. The 2 point calibration is possible using the sequence BUFR\_1-BUFR\_2 or BUFR\_2-BUFR\_3 or even BUFR\_1-BUFR\_3. The 3 point calibration is only possible using the exactly described sequence BUFR\_1-BUFR\_2-BUFR\_3.

		Allowed pH calibration sequences		
r of ion s	1	BUFR_1	BUFR_2	BUFR_3
nbel brat oint	2	BUFR_1 - BUFR_2	BUFR_2 - BUFR_3	BUFR_1 - BUFR_3
Nur cali	3	BUFR_1 - BUFR_2 - BUFR_3		

- A buffer can be disabled in the option selection phase (MENU mode) by choosing the **NIL** value. In this case, the disabled buffer is excluded from the sequence and will not be proposed.
- If no other buffer is expected after the first, as the BUFR\_2 and BUFR\_3 have been set to NIL, the calibration mode is exited by pressing MENU.
- In any case where no operation is performed, the calibration mode automatically stops after 60 seconds.
- If the calibration is rejected by the instrument because it is considered to be excessively corrupted, the **CAL ERR** message will appear, followed by a long beep. The instrument remains in calibration mode and maintains the previous calibration values: at this point, if the calibration is interrupted using ESC, the instrument signals the anomaly through the CAL message blinking.

# Temperature characteristics of Delta OHM pH standard solutions

The 13 standard solutions reported in the table on page 18 are memorized in the instruments with relevant variations according to temperature: The characteristics of the three Delta OHM standard buffers at 6.86 pH, 4.01 pH and 9.18 pH (@25 °C) are reported below.

#### 6.86 pH @ 25 °C

°C	рН	°C	рН
0	6.98	50	6.83
5	6.95	55	6.83
10	6.92	60	6.84
15	6.90	65	6.85
20	6.88	70	6.85
25	6.86	75	6.86
30	6.85	80	6.86
35	6.84	85	6.87
40	6.84	90	6.88
45	6.83	95	6.89

#### 4.01 pH @ 25 °C

°C	Hq	°C	рН
0	4.01	50	4.06
5	4.00	55	4.07
10	4.00	60	4.09
15	4.00	65	4.10
20	4.00	70	4.13
25	4.01	75	4.14
30	4.01	80	4.16
35	4.02	85	4.18
40	4.03	90	4.20
45	4.05	95	4.23

# 9.18 pH @ 25°C

°C	рН	°C	рН
0	9.46	50	9.01
5	9.39	55	8.99
10	9.33	60	8.97
15	9.28	65	8.94
20	9.22	70	8.92
25	9.18	75	8.90
30	9.14	80	8.88
35	9.10	85	8.86
40	9.07	90	8.85
45	9.04	95	8.83

#### **CONDUCTIVITY MEASUREMENT**

The instruments work with conductivity/temperature combined probes, 4-ring and 2-ring, only conductivity probes, or temperature probes. Pt100 or Pt1000 probes may be used for measuring temperature, which is used for the automatic compensation of the conductivity.

The instrument obtains the following from the measurement of conductivity:

- the liquid resistivity measurement ( $\Omega$ , k $\Omega$ , M $\Omega$ ),
- the concentration of total dissolved solids (TDS) according to the X/TDS conversion factor, which can be modified using the menu,
- the salinity (NaCl quantity in the solution, expressed in g/l).

The conductivity, resistivity, TDS or salinity indication is displayed in the main line by pressing **X-\Omega-TDS/ESC**; the secondary line shows the temperature.

The conductivity probes must be periodically calibrated. To facilitate this operation, four automatic calibration solutions recognized by the instrument are provided:

- 0.001 Molar KCI solution (147µS/cm @25°C),
- 0.01 Molar KCl solution (1413µS/cm @25°C),
- 0.1 Molar KCl solution (12880μS/cm @25°C),
- 1 Molar KCl solution (111800µS/cm @25°C),


#### There is no user temperature calibration.

The probes are detected during turn on, and this cannot be performed when the instrument is already on, therefore if a probe is connected and the instrument is on, it must be turned off and back on again.

#### Standard probe

The standard 4-ring conductivity/temperature combined probe has code **SP06T**.

The cell measurement zone is delimited by a bell in PBT. A positioning key, present in the probe's end part, orients the bell correctly when the probe is introduced. For cleaning, simply pull the bell along the probe's axis without rotating it. It is not possible to perform measurements without this bell.

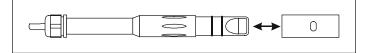

This probe's temperature measuring range is 0...+90 °C.

#### 4-ring or 2-ring probes

The HD2156.1 and HD2156.2 conductivity meters use 4-ring or 2-ring probes for conductivity measurement.

The 4-ring probes are preferred to measure high conductivity solutions, either over an extended range or in presence of pollutants. The 2-ring probes operate in a shorter measurement range but with an accuracy comparable with the 4-ring probes.

The probes can be in glass or plastic: the first can work in presence of aggressive pollutants, the latter are more resistant to collisions, and so more suitable for industrial use.

#### Cells with temperature sensor

The probes fitted with built-in Pt100 or Pt1000 temperature sensor simultaneously measure the conductivity and the temperature: that allows automatic correction of the effect of the latter on the solution conductivity.


# Choosing the cell constant

The constant is a piece of information which characterizes the cell. It depends on its geometry and is expressed in  $cm^{-1}$ . There is no cell capable of measuring the entire conductivity scale accurately enough. Consequently, cells with different constants are used allowing exact measurements on different scales. The cell with constant K = 1 cm<sup>-1</sup> allows measurements from low conductivity up to relatively high conductivity.

The theoretical measurement cell is made of two 1 cm<sup>2</sup> metallic plates separated one from the other by 1 cm. This type of cell has a cell constant Kcell of 1 cm<sup>-1</sup>. In essence, the number, form, material and dimensions of the plates are very different from model to model, from manufacturer to manufacturer.

The low constant K probes are preferably used for low conductivity values, the high constant ones for high values.

The indicative measurement range is reported in the following diagram:

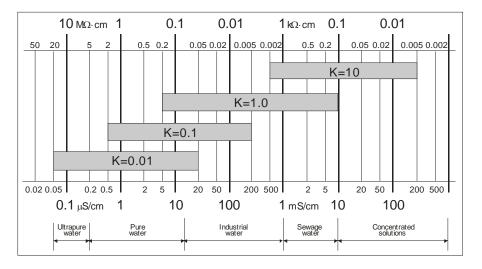

# AUTOMATIC OR MANUAL CONDUCTIVITY COMPENSATION


The conductivity measurement refers to a standard temperature, called reference temperature T\_REF\_COND: that is, the instrument proposes the conductivity you would get if the temperature were T\_REF\_COND. This temperature could be 20 °C or 25 °C according to the setting in the T\_REF\_COND menu item.

The conductivity increase per each grade of temperature variation is a characteristic of the solution and is indicated by the term "temperature coefficient  $\alpha_T$ ": admissible values from 0.00 to 4.00%/°C, default value 2.00%/°C.

When a combined probe with temperature sensor is present, the instrument automatically applies the temperature compensation function, and proposes the measurement using the reference temperature T\_REF according to the coefficient  $\alpha_T$  on the display.

In absence of the temperature probe, the lower display shows the manually set compensation temperature (default=25°C).

To point this condition out, the °C or °F symbol blinks intermittently near the temperature value. On the main display an "m" (manual) is turned on near the battery symbol (if on). The **MT** indication is printed on the print-outs. On the other hand, if the temperature probe is present, the **AT** symbol appears.

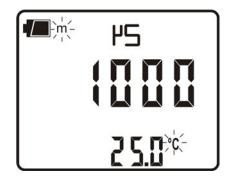

To manually change the compensation temperature, press the °C/°F key once: the indicated temperature value starts blinking. Select the desired temperature value by using the arrows and confirm with ENTER. The display stops blinking, and the temperature displayed is used for compensation.

To change the unit of measurement between °C and °F, press the °C/°F key twice.

#### CALIBRATION OF CONDUCTIVITY

The probe calibration can be carried out on one, two or three points using the standard solutions automatically detected by the instrument (automatic calibration) or other solutions with known value (manual calibration).

The new calibration deletes the data of the previous calibration. For example, if the instrument is calibrated in 4 points and a new 1 point calibration is performed, the instrument will be calibrated in 1 point only.

The CAL symbol will blink when the cell constant is modified using the menu (please see the description of the K\_CELL menu item on page 10).

# Automatic calibration of conductivity using memorized standard solutions

The instrument can recognize four standard calibration solutions:

- 0.001 Molar KCl solution (147µS/cm @25°C),
- 0.01 Molar KCl solution (1413µS/cm @25°C),
- 0.1 Molar KCl solution (12880µS/cm @25°C),
- 1 Molar KCl solution (111800µS/cm @25°C),

Using one of these solutions, the calibration is automatic; the procedure can be repeated with the remaining standard solutions.

The manual calibration is possible with a different conductivity solution from that used in the automatic calibration.

# The solution temperature for the automatic calibration must be between 15°C and 35°C: if the solution temperature is under 15°C or over 35°C, the calibration is rejected: CAL ERR indication appears.

- 1. Turn the instrument on with the ON/OFF key.
- 2. Select the conductivity measurement by pressing "X- $\Omega$ -TDS/ESC".
- 3. Set the probe's cell constant by selecting it from the admitted values: 0.01, 0.1, 0.7, 1.0 or 10.0
- 4. Dip the conductivity meter cell in the calibration solution until the electrodes are covered with liquid.
- 5. Stir the probe lightly to remove any possible air inside the measurement cell.

- 6. If the conductivity probe is not fitted with temperature sensor, press °C/°F and, using the arrows, enter the sample solution temperature value manually (manual setting of temperature). Confirm by pressing ENTER.
- 7. Press the CAL key. The unit of measurement ( $\mu$ S/cm or mS/cm) appears on the comment line. The central line shows the solution conductivity value at the measured temperature, or if the probe is not present, at the manually-set temperature. In the lower line, the closest temperature compensated standard buffer value.

#### If the measurement is in TDS, resistivity or salinity, by pressing CAL, the instrument goes automatically into conductivity calibration mode.

- 8. Press ENTER to confirm the displayed value. The cell constant nominal value (KCELL) and the set temperature coefficient  $\alpha_T$  are displayed. Pressed repeatedly the ENTER key allows the calibration on the point to be repeated, for example, in order to obtain a more stable value.
- 9. To end the probe calibration, press X- $\Omega$ -TDS/ESC.
- 10. Rinse the probe with water. If you are then going to perform low conductivity measurements, we recommend rinsing the probe using distilled or bidistilled water.

The instrument is calibrated and ready for use.

#### Manual calibration of conductivity using non memorized standard solutions

Manual calibration is possible at any calibration solution and temperature if it is within the instrument measurement limits and provided that you know the solution's conductivity at the temperature at which the calibration is performed. Proceed as follows:

- 1. Turn the instrument on with the ON/OFF key.
- 2. Select the conductivity measurement by pressing "X- $\Omega$ -TDS/ESC".
- 3. Set the probe's cell constant by selecting it from the admitted values: 0.01, 0.1, 0.7, 1.0 or 10.0
- 4. Dip the conductivity meter cell into a known conductivity solution until the electrodes are covered with liquid.
- 5. Stir the probe lightly to remove any possible air inside the measurement cell.
- 6. Press MENU, and then FUNC/ENTER until the item ALPH appears. The temperature coefficient  $\alpha_T$  is displayed. Note down the value displayed as it must be set again at the end of the procedure. Set the value to **0.00**. This excludes the temperature compensation during the conductivity measurement.
- 7. Measure the temperature by pressing °C/°F. According to the temperature detected, determine the calibration solution conductivity using the table specifying the conductivity according to temperature.
- 8. Select the conductivity measurement by pressing  $X-\Omega$ -TDS/ESC.
- 9. Press the CAL key. The CAL symbol is turned on. The unit of measurement (μS/cm or mS/cm) appears on the comment line. If the calibration solution conductivity is sufficiently close (-30% to +50%) to one of the standard solutions, the secondary line displays the value. Otherwise it displays the calculated value according to current settings. In the central line the solution conductivity value is indicated according to the cell constant current settings.
- 10. Use the arrows to select the conductivity value determined at point 4 and confirm using ENTER. If the ERR indication appears, see the note below.

- 11. The cell constant nominal value (K\_CELL) and the temperature coefficient  $\alpha_T$  set to 0 are displayed. Pressed repeatedly the ENTER key allows the calibration on the point to be repeated, for example, in order to obtain a more stable value.
- 12. To end the probe calibration, press X- $\Omega$ -TDS/ESC.
- 13. Go back to the MENU and select ALPH: re-enter the temperature coefficient as it was before the calibration.
- 14. Rinse the probe with water. If you are then going to perform low conductivity measurements, we recommend rinsing the probe using distilled or bidistilled water.

The instrument is now calibrated and ready for use.

#### NOTES:

- Without having pressed ENTER at all, the calibration is interrupted by pressing X-Ω-TDS/ESC; the previous values will continue to be used.
- Upon confirming the calibration by using ENTER, the instrument checks that the correction to the conductivity does not exceed the 70% or 150% limits of the theoretical value. If the calibration is rejected because it is considered to be excessively corrupted, the **CAL ERR** message will appear, followed by a long beep. The instrument remains in calibration mode and maintains the previous calibration values: at this point, if the calibration is interrupted using ESC, the instrument signals the anomaly through the CAL message blinking.
- The most frequent causes of error are due to the probe malfunctioning (deposits, dirt...) or to the standard solutions deterioration (bad preservation conditions, alteration due to pollution with different conductivity solutions,...).
- If the measurement is in TDS, resistivity or salinity, on the pressing of CAL, the instrument goes automatically in conductivity calibration.

# Table of standard solution at 147 $\mu\text{S/cm},$ 1413 $\mu\text{S/cm},$ 12.88 mS/cm and 111800 mS/cm

The table reports the standard solutions automatically detected by the instrument according to temperature.

°C	µS/cm	µS∕cm	mS/cm	mS/cm
15.0	121	1147	10.48	92.5
16.0	124	1173	10.72	94.4
17.0	126	1199	10.95	96.3
18.0	128	1225	11.19	98.2
19.0	130	1251	11.43	100.1
20.0	133	1278	11.67	102.1
21.0	136	1305	11.91	104.0
22.0	138	1332	12.15	105.9
23.0	141	1359	12.39	107.9
24.0	144	1386	12.64	109.8
25.0	147	1413	12.88	111.8

°C	µS/cm	µS∕cm	mS/cm	mS/cm
26.0	150	1440	13.13	113.8
27.0	153	1467	13.37	115.7
28.0	157	1494	13.62	117.7
29.0	161	1521	13.87	119.8
30.0	164	1548	14.12	121.9
31.0	168	1581	14.37	124.0
32.0	172	1609	14.62	126.1
33.0	177	1638	14.88	128.3
34.0	181	1667	15.13	130.5
35.0	186	1696	15.39	132.8

#### DIRECT INPUT INTO Pt100 AND Pt1000 TEMPERATURE PROBES

The instrument accepts the input of Platinum temperature probes with resistances of  $100\Omega$  and  $1000\Omega$ .

The Pt100 probes have 4-wire connection; the Pt1000 probes have 2 or 4-wire connection; the excitation current is chosen to minimize the sensor self-heating effects.

The temperature probes are automatically detected by the instrument (please see the description of the Probe Type menu on page 13).

The °C or °F unit of measurement can be chosen for display, printing, and logging using the °C/°F key.

#### How to measure

The temperature measurement by **immersion** is carried out by inserting the probe in the liquid for at least 60 mm; the sensor is housed in the end part of the probe.

In the temperature measurement by **penetration** the probe tip must be inserted to a depth of at least 60mm, the sensor is housed in the end part of the probe. When measuring the temperature on frozen blocks it is convenient to use a mechanical tool to bore a cavity in which to insert the tip probe.

In order to perform a correct **contact** measurement, the measurement surface must be even and smooth, and the probe must be perpendicular to the measurement plane.

So as to obtain the correct measurement, the insertion of a drop of oil or heatconductive paste is useful (do not use water or solvents). This method also improves the response time.

Instructions to connect the TP47 module for conductivity/temperature combined probes, 4-wire Pt100 probes, 2 or 4-wire Pt1000 probes

All Delta Ohm probes are provided with a connector. The HD2156.1 and HD2156.2 instruments also work with third-party conductivity/temperature combined probes, direct 4-wire Pt100 probes, 2 or 4-wire Pt1000 probes. For the instrument connection is available the TP47 connector to which the probe's wires should be welded.




The instructions to connect the probe to the module TP47 are provided below.

The module is supplied complete with fairlead and gasket for 5 mm maximum diameter cables.

Do the following to open the module and connect a probe:

Unscrew the fairlead and extract the gasket, remove the label using a cutter, unscrew the ring on the opposite side as illustrated in the figure:

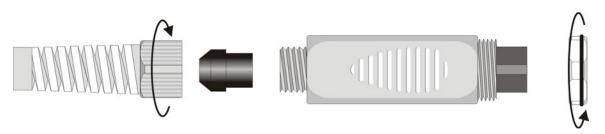

Open the two module shells: the printed circuit to which the temperature probe must be connected is housed inside. The wires coming from the conductivity cell are welded directly onto the pins 1 - 2 - 4 - 5 of the connector.

On the left there are the 1...4 points on which the Pt100 or Pt1000 sensor wires must be welded. The JP1...JP4 jumpers are in the center of the board. These must be left open:

Not Used

Before welding, pass the probe cable through the fairlead and gasket. Weld the temperature sensor wires as shown in the table:

Sensor	TP47 board connection	Jumper
Pt100 4 wires	Pt100 4 wires 4 JP4 4 wires 2 JP3 1 JP1	None
Pt1000 2 wires	Pt1000 2 wires Pt1000 2 wires	JP2
Pt1000 4 wires	Pt1000 4 wires 2 1 JP2 1 JP2	JP2

The wires coming from the **conductivity probe** are welded directly on the DIN45326 connector as reported in the following table:

Sensor	Direct connection to the connector DIN45326	Jumper
2-ring conductivity probe	Internal view of the connector	Jumper between pins 1 and 4 Jumper between pins 2 and 5
4- ring conductivity probe	Internal view of the connector	None


The **Pt100 and Pt1000 sensors** can be welded directly on the DIN45326 connector's pins, as reported in the following table:

Sensor	Direct connection to the connector DIN45326	Jumper
Pt100 4 wires	Pt100 Pt	None
Pt1000 2 wires	Pt1000 Pt1000	Jumper between pins 6 and 7 Jumper between pins 3 and 8

The wires coming from the **conductivity/Pt100 temperature combined probe** are welded directly on the connector as reported in the following table.

Sensor	Direct connection to the connector DIN45326	Jumper
2-ring conductivity probe Pt100 temperature sensor	Pt100 Pt10 Pt1	Jumper between pins 1 and 4 Jumper between pins 2 and 5
4-ring conductivity probe Pt100 temperature sensor	Pt100 Pt	None

Ensure the welds are clean and perfect. Once the welding operation is complete, close the two shells, insert the gasket in the module and screw the fairlead. At the other end of the module, enter the ring with the O-Ring as indicated in the picture.

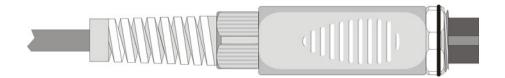

Make sure the cable is not twisted while you are screwing the fairlead. Now the probe is ready.

# WARNINGS AND OPERATING INSTRUCTIONS

- 1. Do not expose the probes to gases or liquids that could corrode the material of the sensor or the probe itself. Clean the probe carefully after each measurement.
- 2. Do not bend the probe connectors or force them upward or downward.
- 3. Do not bend or force the contacts when inserting the probe connector into the instrument.
- 4. Do not bend, deform or drop the probes, as this could cause irreparable damage.
- 5. Always select the most suitable probe for your application.
- 6. Do not use probes in presence of corrosive gases or liquids. The sensor container is made of AISI 316 stainless steel, while the contact probe container is made from AISI 316 stainless steel plus silver. Avoid contact between the probe surface and any sticky surface or substance that could corrode or damage it.
- 7. Above 400°C and below –40°C, avoid violent blows or thermal shocks to Platinum temperature probes as this could cause irreparable damage.
- 8. To obtain reliable measurements, temperature variations that are too rapid must be avoided.
- 9. Temperature probes for surface measurements (contact probes) must be held perpendicular against the surface. Apply oil or heat-conductive paste between the surface and the probe in order to improve contact and reduce reading time. Whatever you do, do not use water or solvent for this purpose. A contact measurement is always very hard to perform. It has high levels of uncertainty and depends on the ability of the operator.
- 10. Temperature measurements on non-metal surfaces usually require a great deal of time due to the low heat conductivity of non-metal materials.
- 11. Probes are not insulated from their external casing; be very careful not to come into contact with live parts (above 48V). This could be extremely dangerous for the in-



strument as well as for the operator, who could be electrocuted.

12. Avoid taking measurements in presence of high frequency sources, microwave ovens or large magnetic fields; results may not be very reliable.

- 13. Clean the probe carefully after use.
- 14. The instrument is water resistant and IP66, but is not watertight and therefore should not be immersed in water without closing the free connectors using caps. **The probe connectors must be fitted with sealing gaskets.** Should the instrument fall into the water, check for any water infiltration. Gently handle the instrument in such a way as to prevent any water infiltration from the connectors' side.

#### Maintenance

The instrument case is made of ABS plastic and the protections are rubber: do not use any incompatible solvent for cleaning.

Do not use any abrasive products for cleaning.

#### NOTES ON THE USE OF THE **pH** ELECTRODES

The average life of a pH electrode is about 1 year, depending on the use and maintenance.

The electrodes used at high temperatures or in strong alkaline environments normally have a shorter life.

The new electrodes should be conditioned for half a day immersed in a 6.86pH or 4pH buffer.

Calibrate the electrode using buffer solutions close to the values to be measured. A new electrode should always be calibrated at neutral pH (6.86pH) as first point, and at least in a second point.

Following are listed some problems that frequently occur and the possible solutions.

Wrong pH measure. Perform the following checks:

Check the diaphragm is not blocked and if necessary use a cleaning solution.

Check the reference system is not polluted and, if the electrode is refillable, replace the electrolyte.

Check that the tip of the electrode has no air bubbles inside and that the electrode is sufficiently immersed.

Even residues of dirt deposited on the membrane can alter the measure: use a solution for protein cleaning.

**Slow response and wrong measures**. A possible cause can be the aging or the erosion of the membrane, or a short-circuit in the connector.

**Storage.** Store the electrode immersed in a maintenance solution specific for pH electrodes.

#### NOTES ON CONDUCTIVITY PROBES

The service life of a cell can be unlimited, provided that the necessary maintenance is performed and that it does not break. Some of the most frequent problems and their possible solutions are reported below.

**Measurement of conductivity different from the expected value**. Check that the cell used is suitable for the measurement range. Check that the cell is not dirty, that there are no air bubbles inside it. Calibrate again using the appropriate standard.

**Slow response or instability**. Check that the cell is not dirty, that there are no traces of oil or air bubbles inside it. If you work with a black Platinum cell, new regeneration of the electrode could be necessary.

**Cell constant value not accepted**. Check that the standard solutions are in good condition, that the probe's constant cell value coincides with that selected in the instrument and that the solution temperature is within the range 15...35 °C.

# **INSTRUMENT SIGNALS AND FAULTS**

The following table lists all error indications and information displayed by the instrument and supplied to the user in different operating situations:


Display indications	Explanation	
	This appears if the conductivity/temperature probe is measuring a value exceeding the set measuring range.	
ERR	This message appears if the pH measurement exceeds the -2.00pH19.99pH limits, if the mV measurement exceeds the $\pm 2.4V$ limits, if the temperature probe, already detected by the instrument, is disconnected.	
CAL	This appears when, during calibration, the read value exceeds the limits of $-30\%$ or $+50\%$ of the standard solution value compensated for temperature or the solution temperature is less than $15^{\circ}$ C or higher than $35^{\circ}$ C.	
ERR	In pH when the offset and/or slope limits exceed the following values:  Offset  > 20mV Slope < 50mV/pH or Slope > 63mV/pH.	
LOG MEM FULL	Memory full; the instrument cannot store further data, the memory space is full.	
CAL blinking	No calibration has yet been effected, or the cell constant value has been modified in the menu (please see the description of the K_CELL menu item on page 12).	
	In pH the calibration was not completed correctly.	
m	No probe with temperature sensor connected. The letter "m" indicates that the displayed temperature was entered manually.	
OVER	Measurement overflow: this appears when the probe measures a value exceeding the measurement range or the mV measurement is included in the +2.0+2.4V range.	
UNDR	Measurement overflow: this appears when the mV measurement is included in the -2.42.0V range.	
SYS ERR #	Instrument management program error. Contact the instrument's supplier and communicate the numeric code # reported by the display.	
CAL LOST	Program error: it appears after turning on for a few seconds. Contact the instrument's supplier.	
BATT TOO LOW CHNG NOW	Insufficient battery charge appearing on turning on. The instrument issues a long beep and turns off. Replace the batteries.	

The following table reports the indications provided by the instrument as they appear on the display, together with their description.

Display indication	Explanation		
ALPH	temperature coefficient $\alpha_T$		
AUTO-HOLD	automatic function holding the displayed meas- urement		
BATT TOO LOW - CHNG NOW	battery discharged - replace it immediately		
BAUDRATE	baud rate value		
BUFR_1_pH	first pH buffer value		
BUFR_2_pH	second pH buffer value		
BUFR_3_pH	third pH buffer value		
CAL ERR	calibration error		
COMM STOP	printing complete		
COMM STRT	printing started		
DAY_	day		
DUMP_END	data transfer complete		
DUMP_In_PROG	data transfer in progress		
FUNC CLR	max, min and average values clearing		
FUNC CLRD	max, min and average values cleared		
HOUR	hour		
KCEL_COND	conductivity cell constant K		
КОНМ	measurement of resistivity in $k\Omega$		
LAST CAL COND m/d h/m	last conductivity calibration date month/day hour/minutes		
LAST CAL pH m/d h/m	last pH calibration date month/day hour/minutes		
LOG IN PROG	logging in progress		
LOG MEM FULL	memory full		
LOG_CLRD	memory data cleared		
LOG_DUMP_OR_ERAS	transfer or erase data		
LOG_STOP	logging complete		
LOG_STRT	logging started		
MIN >>> USE_UNIT_TO_ZERO SEC	minutes >>> use the UNIT key to reset the sec- onds		
MNTH	month		
МОНМ	measurement of resistivity in $M\Omega$		
NaCl	measurement of salinity in g/l		
NIL	disable current pH buffer		
OFFS	pH measurement offset		
ОНМ	measurement of resistivity in $\Omega$		
OVER	maximum expected range limit exceeded		
PLS_EXIT >>> FUNC RES_FOR_FACT	please exit using ESC >>> function reserved to		
ONLY	factory calibration		
PRBE_TYPE	type of probe connected		
PRNT AND LOG INTV	printing and logging intervals		
PRNT INTV	printing interval		
PROB COMM LOST	lost communication with probe		
PROB ERR	error - unexpected probe		
RCD MODE	record function operating mode (max, min, avg)		
REF_TEMP COND	conductivity reference temperature		
SET USER BUFR pH	user defined pH buffer value		
SLP_MODE_LOG	turning off during recording mode		
SLPE%	gain in % during pH calibration		
SMPL ID REL=RSET SER=PRINT	identifier of the sample - REL=reset -		

Display indication	Explanation
	SERIAL=print heading
SYS ERR #	program error number #
TDS	total dissolved solids
UNDR	minimum expected range limit exceeded
YEAR	year

# LOW BATTERY WARNING AND BATTERY REPLACEMENT

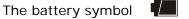

on the display constantly shows the battery charge status. To the extent that batteries have discharged, the symbol "empties". When the charge decreases still further it starts blinking...



In this case, batteries should be replaced as soon as possible.

If you continue to use it, the instrument can no longer ensure correct measurement. The memory data are maintained.

If the battery charge level is insufficient, the following message appears when you turn the instrument on:

#### BATT TOO LOW CHNG NOW

The instrument issues a long beep and turns off. In this case, replace the batteries in order to turn the instrument back on.

In order to avoid data loss, the logging session is ended, if the HD2156.2 is logging and battery voltage falls below the minimum operating level.

The battery symbol turns off when the external power supply is connected.

To replace the batteries, switch the instrument off and unscrew the battery cover counter clockwise. After replacing the batteries (4 1.5V alkaline batteries - type AA) screw the cover on clockwise.



After replacing the batteries, the date, time, baud rate, type of probe, printing interval, logging parameters must be set again: in order to simplify the operation, on insertion of the new batteries the instrument turns on automatically and requests these parameters in sequence. To go to the next item press ENTER; to return to measurement mode, press MENU.

#### MALFUNCTIONING UPON TURNING ON AFTER BATTERY REPLACEMENT

After replacing the batteries, the instrument may not restart correctly; in this case, repeat the operation. After disconnecting the batteries, wait a few minutes in order to allow circuit condensers to discharge completely; then reinsert the batteries.

#### WARNING ABOUT BATTERY USE

- Batteries should be removed when the instrument is not used for an extended time.
- Flat batteries must be replaced immediately.
- Avoid batteries leaking.
- Always use good quality leakproof alkaline batteries. Sometimes on the market, it is possible to find new batteries with an insufficient charge capacity.

# **INSTRUMENT STORAGE**

Instrument storage conditions:

- Temperature: -25...+65°C.
- Humidity: less than 90%RH without condensation.
- Do not store the instrument in places where:

Humidity is high. The instrument may be exposed to direct sunlight. The instrument may be exposed to a source of high temperature. The instrument may be exposed to strong vibrations. The instrument may be exposed to steam, salt or any corrosive gas.

# SERIAL INTERFACE AND USB

The HD2156.1 and HD2156.2 instruments are fitted with an electrically isolated RS-232C serial interface; the HD2156.2 also has an USB 2.0 interface.

The following serial cables can be used:

- HD2110CSNM: serial connection cable with 8-pole MiniDin connector on one end and 9-pole Sub D connector on the other end;
- C.206: serial connection cable with 8-pole MiniDin connector on one end and USB type A connector on the other end. With integrated RS232/USB converter;
- CP23: connection cable with Mini-USB type B connector on one end and USB type A connector on the other end (only for HD2156.2).

The connection via the C.206 cable requires the previous installation of the cable USB drivers. Install the drivers before connecting the C.206 cable to the PC.

The connection via the CP23 cable does not require the installation of USB drivers: when connecting the instrument to the PC, the Windows® operating system automatically recognizes the device as an HID device (Human Interface Device) and uses the drivers already included in the operating system.

	Cable	Instrument port	PC port	Installation of USB drivers
	HD2110CSNM	RS232 (MiniDin)	RS232 (9-pole SubD)	No
Ī	C.206	RS232 (MiniDin)	USB	Yes
	CP23	USB (Mini-USB)	USB	No

The instrument standard serial transmission parameters are:

Baud rate 38400 baud

8

1

- Parity None
- N. bit
- Stop bit
- Protocol Xon/Xoff

It is possible to change the RS232C serial port baud rate by setting the "Baudrate" parameter in the menu (please see page 13). The possible values are: 38400, 19200, 9600, 4800, 2400, 1200. The other transmission parameters are fixed.

The USB 2.0 connection does not require the setting of parameters.

The instruments are provided with a complete set of commands and data queries to be sent via the PC. The serial commands work with a standard serial communication program (e.g. Hyperterminal) only through the RS232 serial port of the instrument, using the cable HD2110CSNM or the cable C.206.

All the commands transferred to the instrument must have the following structure: **XYcr** where: **XY** is the command code and **cr** is the Carriage Return (ASCII 0D)

Com- mand	Response	Description
PO	&	Ping (locks the instrument keyboard for 70 seconds)
P1	&	Unlocks the instrument keyboard
S0		Captured measurements (24 characters) AT = automat- ic temperature compensation, temperature
G0 Model HD2156 -2		Instrument model
G1	M=pH/conductivity meter	Model description

Com- mand	Response	Description	
G2	SN=12345678	Instrument serial number	
G3	Firm.Ver.=01-01	Firmware version	
G4	Firm.Date=2004/06/15	Firmware date	
G5	cal 0000/00/00 00:00:00	Calibration date and time	
G6	Probe A= Cond.+Pt1000	Type of probe connected to input	
GB	User ID=000000000000000000000000000000000000	User code (set with T2xxxxxxxxxxxxxxxxx)	
GC		Print instrument's heading	
КО		Stop printing data	
K1	PRINTOUT IMMEDIATE MODE	Immediate printing of data	
К4	&	Start logging data	
К5	&	Stop logging data	
КР	&	Auto-power-off function=ENABLE	
KQ	&	Auto-power-off function=DISABLE	
LD	PRINTOUT OF LOG	Print data logged in flash	
LE	&	Erase data in flash memory	
LN	&2000	Number of free pages in the flash memory	
LUAn	&	Sets the unit of measurement for the temperature n=0 > °C n=1 > °F	
LUBn	&	Sets the unit of measurement for the conductivity $n=0 > \mu S/cm$ $n=1 > \Omega$ n=2 > TDS n=3 > NaCl	
LUCn	&	Sets the unit of measurement for the pH $n=0 > pH$ n=1 > mV	
RA	& #	Reading of LOG/PRINT interval set	
RP	& 700	Battery level (Resolut. 0.01V)	
RUA	U= °C	Temperature measurement unit	
RUB	U= uS	Conductivity measurement unit	
RUC	U= pH	pH measurement unit	
WA#	&	Setting LOG/PRINT interval. # is a hexadecimal number 0D that represents the po- sition of the interval in the list 0, 1, 5, 10,, 3600 sec- onds.	
WC0	&	Setting SELF off	
WC1	&	Setting SELF on	

Command characters are exclusively upper case characters. Once a correct command is entered, the instrument responds with "&"; when any wrong combination of characters is entered, the instrument responds with "?". The instrument response strings end with the sending of the CR command (Carriage Return). The instrument does not send the LF command (Line Feed).

Before sending commands to the instrument via the serial port, locking the keyboard to avoid functioning conflicts is recommended: use the P0 command. When complete, restore the keyboard with the P1 command.

# STORING AND TRANSFERRING DATA TO A PERSONAL COMPUTER

The HD2156.1 and HD2156.2 instruments can be connected to a personal computer via an RS232C serial port or USB 2.0 port, and exchange data and information through the DeltaLog9 software running in a Windows operating environment. Both models can send in real time input measured values directly to a PC, through the PRINT function; the HD2156.2 can also store the values measured by using the *Logging* function (LOG key) in its internal memory. If necessary, the data stored in the memory can be transferred to a PC later.

# THE *LOGGING* FUNCTION - ONLY FOR HD2156.2

The *Logging* function allows the recording up to 20,000 sets of three measurements [T-X-pH] registered by the probes connected to the inputs. Logging always includes **three** parameters. Each set of three data is composed of: temperature in °C or °F, conductivity or resistivity or TDS or NaCl, pH or mV.

The logged parameters are selected using the "°C/°F", "pH/mV-▼" and "X-Ω-TDS/ESC" keys.

The time interval between two consecutive measurements can be set from 1 second to 1 hour. The logging starts by pressing the LOG key and ends by pressing the same key again: the data memorized in this way form a continuous block of samples.

See the description of the menu items on page 10.

If the automatic turning off option between two recordings (MENU >> **Sleep\_Mode\_LOG**) is enabled, upon pressing the LOG key the instrument logs the first data and turns off (the LOG ON indication periodically appears on the display to indicate that the Logging function is active). 15 seconds before the next logging instant, it turns on again to capture the new sample, and then turns off.

The data stored in the memory can be transferred to a PC using the DUMP LOG command: MENU >> LOG. During data transfer the display shows the message DUMP; to stop the data transfer press ESC on the instrument or on the PC.

## CLEARING THE MEMORY - ONLY FOR HD2156.2

To clear the memory, use the Erase Log function (MENU >> SERIAL/Erase Log).

The instrument starts clearing the internal memory; at the end of the operation, it goes back to normal display.

NOTES:

- Data transfer does not cause the memory to be erased; the operation can be repeated as many times as required.
- The stored data remain in the memory independently of battery charge conditions.
- In order to print the data to a parallel interface printer, you must use a parallel-serial adaptor (not supplied).
- The direct connection between instrument and printer via a USB connector does not work.
- Some keys are disabled during logging. The following keys work: ON/OFF, FUNC (Max-Min-Avg) and SERIAL.
- The recording started with the display in Max-Min-Avg mode proceeds normally with the actual measured values. Only the display shows respectively the Max, Min or Avg values.
- The logging is disabled, if the Auto-HOLD function is enabled.
- It is possible to activate both the logging (LOG) and direct transfer (PRINT) functions at the same time.

# THE **PRINT** FUNCTION

The PRINT function sends the measurements taken in real time by the instrument inputs directly to a PC or a printer. Print data units of measurements are the same as those used on the display. Printing always includes **three** parameters. Each set of three data is composed of: temperature in °C or °F, conductivity or resistivity or TDS or NaCl, pH or mV. The printed parameters are selected using the "°C/°F", "pH/mV- $\checkmark$ " and "X- $\Omega$ -TDS/ESC" keys.

The function is started by pressing SERIAL. The time interval between two consecutive prints can be set from 1 second to 1 hour (please see the **Print and log interval** menu item on page 10). If the print interval is equal to 0, by pressing PRINT the single data is sent to the connected device. If the print interval is higher than 0, the data transfer continues until the operator stops it by pressing SERIAL again. See the description of the menu items on page 10.

The PRINT function works with a standard serial communication program only through the RS232 serial port of the instrument, using the cable HD2110CSNM or the cable C.206.

Connect the HD40.1 printer using cable HD2110CSNM.

NOTES:

- The print out is formatted across 24 columns.
- Some keys are disabled during serial transmission. The following keys work: ON/OFF, FUNC (Max-Min-Avg) and LOG.
- The FUNC key has no effect on the print-outs, only on the display.
- If the Auto-Hold function is enabled, the print time interval is forced to zero: pressing SERIAL only causes the print to occur when the measurement has stabilized (HOLD symbol still). Later, it is possible to repeat the print at will, but while the HOLD mode is on, the sample identifier number is not increased. This is useful when more labels must be printed with the same identification code without increasing the code each time.
- It is possible to activate both the logging (LOG) and direct transfer (PRINT) functions at the same time.

INTOUT INMEDIATE MODE del HD2156 -2	\ ſ			
ACONDUCTIVITY METER		Date=2	004/01/1	05:22:23
irm. ver. = 2004/09/24		MT		1001
irm. 046678 N=12345678			25.0 1413	L
N=12345670 Cal.=Factory Jser ID=000000000000000000	0		8.26	u S P H
User ID=00000000		Date=20	84/81/0	F.I.
K.Cell 0.905:03:	35	Sample	ID=0000	1 05:22:44
time 2004 0.00		MT	25.0	00
alpha-' a 50			11.49 7.06	mS
TD5 Temp.Ref. = 25.0°C Temp.Ref. = Pt1000		21- 00-		РH
Probe Lar	8 S	are=206	04/01/01 D=00000	05:22:54
0-1 Temp. = 11 05:03	:14 M	T	25.0	3
time 2004/01/01 05.05 0fs=-13.7 mV Slp=100	. Ø Z		412	°C uS
Ofs=-13. (	9:45		7.06	DH
0fs=-13.7 m # Date=2004/01/01 05:2 Date=interval=	sec Da	te=200	4/01/01	05:23:39
Comple		imple I		
MT 20.0 U	s		25.0 50.5	°C
1414 7.36 F	H		7.06	G/L PH
	Dai	te=2004	104.14	05:24:01
	us San	nple ID	-0000005	35:24:01
7.86	PH MT		25.0	°C
25.8	°C	•	19.7	онм
MT 433	uS pH Dat		7.06	PH
7.86	Sam	e=2004/ ple ID=	01/01 0	5:24:39
25.0	°C MT		0000000	
MT 3.47	mS pH			°C G/L
7.93			1 05	PH
> Printout sto	pped			

### CONNECTION TO A PC

# <u>HD2156.1</u>

Connection to the PC with the cable:

- HD2110CSNM: 8-pole MiniDin connector on one end and 9-pole Sub D connector on the other end;
- C.206: 8-pole MiniDin connector on one end and USB type A connector on the other end. With integrated RS232/USB converter (requires the installation of the USB drivers).

## <u>HD2156.2</u>

Connection to the PC with the cable:

- CP23: Mini-USB type B connector on one end and USB type A connector on the other end;
- HD2110CSNM: 8-pole MiniDin connector on one end and 9-pole Sub D connector on the other end;
- C.206: 8-pole MiniDin connector on one end and USB type A connector on the other end. With integrated RS232/USB converter (requires the installation of the USB drivers).

The instruments are supplied with the DeltaLog9 software that manages the connection, data transfer, graphic presentation, and printing operations of the captured or logged measurements.

The DeltaLog9 software is complete with "On-line Help" (also in PDF format) describing its characteristics and functions.

#### CONNECTION TO THE RS232C SERIAL PORT OF THE INSTRUMENT

- 1. The measurement instrument must be switched off.
- 2. Using the Delta Ohm HD2110CSNM or C.206 cable, connect the measurement instrument to the first free RS232C (COM) or USB serial port of the PC.
- 3. Turn on the instrument and set the baud rate to 38400 (MENU >> ENTER until the Baud Rate parameter >> select 38400 using the arrows >> confirm with ENTER). The parameter remains in the memory until replacement of the batteries.
- 4. Start the DeltaLog9 application and press CONNECT. Wait for the connection to occur and follow the indications on the screen. For a description of the DeltaLog9 application, please refer to its on-line Help.

# CONNECTION TO THE USB 2.0 PORT OF THE INSTRUMENT - ONLY FOR HD2156.2

The connection via the CP23 cable does not require the installation of USB drivers: when connecting the instrument to the PC, the Windows® operating system automatically recognizes the device as an HID device (Human Interface Device) and uses the drivers already included in the operating system.

To check if the connection has been successfully completed, double-click on "*Device Manager*" from the Control Panel. The following items should appear:

"Human Interface Device" >> "HID-compliant device"

"Human Interface Device" >> "USB Human Interface Device"

When the USB cable is disconnected, the items disappear and reappear when it is connected again.

# NOTES ABOUT WORKING AND OPERATIVE SAFETY

## Authorized use

The technical specifications as given in chapter "TECHNICAL CHARACTERISTICS" must be observed. Only the operation and running of the measuring instrument according to the instructions given in this operating manual is authorized. Any other use is considered unauthorized.

# General safety instructions

This measuring system is constructed and tested in compliance with the EN 61010-1:2010 safety regulations for electronic measuring instruments. It left the factory in a safe and secure technical condition.

The smooth functioning and operational safety of the measuring system can only be guaranteed if the generally applicable safety measures and the specific safety instructions in this operating manual are followed during operation.

The smooth functioning and operational safety of the instrument can only be guaranteed under the environmental and electrical operating conditions that are in specified in chapter "TECHNICAL CHARACTERISTICS".

Do not use or store the product in places such as listed below:

- Rapid changes in ambient temperature which may cause condensation.
- Corrosive or inflammable gases.
- Direct vibration or shock to the instrument.
- Excessive induction noise, static electricity, magnetic fields or noise.

If the measuring system was transported from a cold environment to a warm environment, the formation of condensate can impair the functioning of the measuring system. In this event, wait until the temperature of the measuring system reaches room temperature before putting the measuring system back into operation.

## Obligations of the purchaser

The purchaser of this measuring system must ensure that the following laws and guidelines are observed when using dangerous substances:

- EEC directives for protective labour legislation
- National protective labour legislation
- Safety regulations

# **INSTRUMENT TECHNICAL CHARACTERISTICS**

# Instrument

Dimensions (Length x Width x Height) Weight Materials Display

Operating conditions Working temperature Storing temperature Working relative humidity Protection degree

#### Power Supply

Batteries Autonomy Power absorbed with instrument off Mains (cod. **SWD10**)

Security of memorized data

# Time

Date and time Accuracy *Measured values storage - model* **HD2156.2** Type Quantity

# Selectable storage interval

# Serial interface RS232C

Type Baud rate Data bit Parity Stop bit Flow Control Serial cable length Selectable print interval

USB interface - model **HD2156.2** Type 185 x 90 x 40 mm
470g (complete with batteries)
ABS, rubber
2x4½ digits plus symbols
Visible area: 52x42mm

-5...50°C -25...65°C 0...90%RH without condensation IP66

4 1.5V type AA batteries 200 hours with 1800mAh alkaline batteries 20μA Output mains adapter 100-240 Vac/12Vdc-1A Unlimited, independent of battery charge conditions

Schedule in real time 1min/month max drift

2000 pages containing 10 samples each 20000 sets of measures composed of [pH or mV], [X or  $\Omega$  or TDS or NaCl], [°C or °F] depending on the configuration.

1, 5, 10, 15, 30 s, 1min, 2min, 5min, 10min, 15min, 20min, 30min and 1hour

RS232C electrically isolated Can be set from 1200 to 38400 baud 8 None 1 Xon/Xoff Max 15m immediate or 1s, 5s, 10s, 15s, 30s, 1min, 2min, 5min, 10min, 15min, 20min, 30min and 1hour

1.1 - 2.0 electrically isolated

Connections	
pH/mV input	Female BNC connector
Conductivity input	8-pole male DIN45326 connector
RS232 serial interface	8-pole MiniDin connector
USB interface (only HD2156.2)	Mini-USB type B connector
Mains adapter (cod. SWD10)	2-pole connector (positive at centre)
Measurement of pH by Instrument	
Measurement range	-2.00+19.99 pH
Resolution	0.01 pH
Accuracy	±0.01 pH ±1digit
Input impedance	>10 <sup>12</sup> Ω
Calibration error @25°C	Offset  > 20 mV
	Slope > 63 mV/pH or Slope < 50 mV/pH
	Sensitivity > 106.5% or Sensitivity < 85%
Automatic/manual temperature compensatio	n -50+150 °C
Measurement of mV by Instrument	
Measurement range	-1999.9+1999.9 mV
Resolution	0.1 mV
Accuracy	$\pm 1$ mV or $\pm 0.2\%$ of the reading (the greater)
Drift after 1 year	0.5 mV/year
Standard solutions automatically detected (@25°C	)
	1.679 pH - 2.000 pH - 4.000 pH - 4.008 pH - 4.010 pH - 6.860 pH - 6.865 pH - 7.000 pH - 7.413 pH - 7.648 pH - 9.180 pH - 9.210 pH - 10.010 pH
Measurement of conductivity by Instrument	
Measurement range (Kcell=0.01) / Resolutio	n 0.000019.999 μS/cm / 0.001 μS/cm
Measurement range (Kcell=0.1) / Resolution	0.0019.99 μS/cm / 0.01 μS/cm
Measurement range (K cell=1) / Resolution	0.0199.9 μS/cm / 0.1 μS/cm
	2001999 μS/cm / 1 μS/cm
	2.0019.99 mS/cm / 0.01 mS/cm
	20.0199.9 mS/cm / 0.1 mS/cm
Measurement range (K cell=10) / Resolution	
Accuracy (conductivity)	±0.5% ±1digit

Measurement of resistivity by Instrument Measurement range (Kcell=0.01) / Resolutior	$1 \ln t_0 1 G\Omega_{cm} / (*)$
Measurement range (Kcell=0.1) / Resolution	1 , ,
Measurement range (Kcell=1) / Resolution	5.0199.9Ω·cm / 0.1Ω·cm
	200999Ω·cm / 1Ω·cm
	1.00k19.99kΩ·cm / 0.01kΩ·cm
	20.0k99.9kΩ·cm / 0.1kΩ·cm
	100k999kΩ·cm / 1kΩ·cm
	110MΩ·cm / 1MΩ·cm
Measurement range (Kcell=10) / Resolution	0.55.0Ω·cm / 0.1Ω·cm
Accuracy (resistivity)	±0.5% ±1digit
Measurement of total dissolved solids (with coefficient Measurement range (Kcell=0.01) / Resolution	0.00019.999 mg/l / 0.005 mg/l
Measurement range (Kcell=0.1) / Resolution Measurement range (K cell=1) / Resolution	
	20.099.9 g/l / 0.1 g/l
Measurement range (K cell=10) / Resolution	<b>c c</b>
Accuracy (total dissolved solids)	±0.5% ±1digit
Measurement of salinity	0
Measurement range / Resolution	0.0001.999 g/l / 1 mg/l 2.0019.99 g/l / 10 mg/l 20.0199.9 g/l / 0.1 g/l
Accuracy (salinity)	±0.5% ±1digit

(\*) The resistivity measurement is obtained from the reciprocal of conductivity measurement. Close to the bottom of the scale, the indication of resistivity appears like reported in the table below.

K cell = $0.01 \text{ cm}^{-1}$		
Conductivity(µS/cm)	Resistivity (MΩ·cm)	
0.001 µS/cm	1000 MΩ·cm	
0.002 µS/cm	500 MΩ·cm	
0.003 µS/cm	333 MΩ·cm	
0.004 µS/cm	250 MΩ·cm	

K cell = $0.1 \text{ cm}^{-1}$		
Conductivity (µS/cm)	Resistivity (MΩ·cm)	
0.01 µS/cm	100 MΩ·cm	
0.02 µS/cm	50 MΩ·cm	
0.03 µS/cm	33 MΩ·cm	
0.04 µS/cm	25 MΩ·cm	
•••		

(\*) The resistivity measurement is obtained from the reciprocal of conductivity measurement. Close to the bottom of the scale, the indication of resistivity appears like reported in the table below.

K cell = $0.01 \text{ cm}^{-1}$		
Conductivity(µS/cm)	Resistivity (MΩ·cm)	
0.001 µS/cm	1000 MΩ·cm	
0.002 µS/cm	500 MΩ·cm	
0.003 µS/cm	333 MΩ·cm	
0.004 µS/cm	250 MΩ·cm	

K cell = 0.1 cm <sup>-1</sup>		
Conductivity (µS/cm)	Resistivity (MΩ·cm)	
0.01 µS/cm	100 MΩ·cm	
0.02 µS/cm	50 MΩ·cm	
0.03 µS/cm	33 MΩ·cm	
0.04 µS/cm	25 MΩ·cm	

Automatic/manual temperature compensation	0100 °C with $\alpha_T$ that can be selected from 0.00 to 4.00%/°C
Reference temperature	20 °C or 25 °C
X/TDS conversion factor	0.40.8
Cell constant K (cm <sup>-1</sup> )	0.01, 0.1, 0.7, 1.0 and 10.0
Standard solutions automatically detected (@25°C	<u>;)</u>
	147 μS/cm
	1413 μS/cm
	12880 μS/cm
	111800 μS/cm
Measurement of temperature by Instrument	
Pt100 measurement range	-50+200 °C
Pt1000 measurement range	-50+200 °C
Resolution	0.1 °C
Accuracy	±0.25 °C
Drift after 1 year	0.1 °C/year

# **TECHNICAL DATA OF PROBES**

# pH electrodes

ORDERING CODE	MEASURING RANGE AND USE	DIMENSIONS
КР30	014 pH / 080 °C / 2bar Glass body - GEL filled 1 ceramic diaphragm Cable L=1 m with BNC Waste water, Drinking water, Water emul- sions, Galvanic, Paints, Varnish, Water sus- pensions, Fruit-juices Titration.	Ø 16 130 Ø 12 BNC

# 2 AND 4-ELECTRODE CONDUCTIVITY PROBES

ORDER CODE	MEASUREMENT RANGE	DIMENSIONS
SP06T	K=0.7 5 μS/cm100 mS/cm 090 °C 4-electrode cell PBT/Platinum	L=1.5m
SPT01G	K=0.1 0.1 μS/cm500 μS/cm 080 °C 2-electrode cell Glass/Platinum	L=1.5m
SPT1G	K=1 10 μS/cm10 mS/cm 080 °C 2-electrode cell Glass/Platinum	L=1.5m
SPT10G	K=10 500 µS/cm200 mS/cm 080 °C 2-electrode cell Glass/Platinum	L=1.5m

# Pt100 4-WIRE PROBES AND Pt1000 2-WIRE PROBES COMPLETE WITH TP47 MODULE

Model	Туре	Application range	Accuracy
TP47.100.0	Pt100 4 wires	-50+250°C	Class A
TP47.1000.0	Pt1000 4 wires	-50+250°C	Class A
TP87.100.0	Pt100 4 wires	-50+200°C	Class A
TP87.1000.0	Pt1000 4 wires	-50+200°C	Class A

Common characteristics

# Resolution

#### 0.01°C nel campo ±199.99°C,

0.1°C altrove

Temperature drift @ 20°C Pt100 0.003%/°C Pt1000 0.005%/°C

# **ORDERING CODES**

- **HD2156.1** Kit including: instrument HD2156.1, 4 x 1.5V alkaline batteries, operating manual, case and DeltaLog9 software (downloadable from Delta OHM website).
- HD2156.2 Kit including: instrument HD2156.2 data logger, 4 x 1.5V alkaline batteries, operating manual, CP23 USB cable, case and DeltaLog9 software (down-loadable from Delta OHM website).

pH/mV electrodes, conductivity probes, temperature probes, standard reference solutions, connection cables for pH electrodes with S7 connector, HD2110CSNM and C.206 cables have to be ordered separately.

- HD2110CSNM Connection cable 8-pole MiniDin Sub D 9-pole female for RS232C.
- C.206 Connection cable 8-pole MiniDin USB type A. With integrated RS232/USB converter.
- **SWD10** Stabilized power supply at 100-240 Vac/12 Vdc-1A mains voltage.
- **HD40.1** The kit includes: 24-column portable thermal printer, serial interface, 57 mm paper width, four NiMH 1.2V rechargeable batteries, SWD10 power supply, instruction manual, 5 thermal paper rolls.
- **BAT-40** Spare battery pack for HD40.1 printer with in-built temperature sensor.
- **RCT** The kit includes 4 thermal paper rolls 57 mm wide and 32 mm in diameter.
- HD22.2 Laboratory electrode holder composed of basis plate with incorporated magnetic stirrer, staff and replaceable electrode holder. Height max. 380 mm. For Ø12 mm electrodes.
- **HD22.3** Laboratory electrode holder with metal basis plate. Flexible electrode holder for free positioning. For Ø12 mm probes.

#### pH ELECTRODES

- **KP30** Combined pH electrode for common use, GEL filled, cable 1m with BNC, Epoxy body
- **CP** Extension cable 1.5 m with BNC connector on one side and S7 on the other side, for electrode without cable with S7 connector.
- **BNC** Female BNC for electrode extension.

#### pH STANDARD SOLUTIONS

**HD8600C** ACCREDIA ISO 17025 certified buffer solutions kit: pH 4.01+ pH 7.00 + pH 10.01. Two 50 ml bottles for each type (6 bottles in total).

#### CONDUCTIVITY PROBES AND COMBINED CONDUCTIVITY AND TEMPERATURE PROBES

- **SP06T** Combined conductivity and temperature 4-electrode cell in Platinum, body in PBT. Cell constant K = 0.7. Measurement range 5  $\mu$ S/cm...100 mS/cm, 0...90 °C.
- **SPT01G** Combined conductivity and temperature 2-electrode Platinum-wire cell, glass body. Cell constant K = 0.1. Measurement range 0.1  $\mu$ S/cm...500  $\mu$ S/cm, 0...80 °C.
- **SPT1G** Combined conductivity and temperature 2-electrode Platinum-wire cell, glass body. Cell constant K = 1. Measurement range 10  $\mu$ S/cm...10 mS/cm, 0...80 °C.

**SPT10G** Combined conductivity and temperature 2-electrode Platinum-wire cell, glass body. Cell constant K = 10. Measurement range 500  $\mu$ S/cm...200 mS/cm, 0...80 °C.

#### STANDARD CONDUCTIVITY CALIBRATION SOLUTIONS

HD8700CACCREDIA ISO 17025 standard solutions kit: 0.001 mol/l (147  $\mu$ S/cm @25<br/>°C) + 0.01 mol/l (1413  $\mu$ S/cm @25 °C) + 0.1 mol/l (12880  $\mu$ S/cm @25 °C).<br/>Two 50 ml bottles for each type (6 bottles in total).

#### TEMPERATURE PROBES COMPLETE WITH TP47 MODULE

- **TP47.100.0** Immersion probe, sensor Pt100 direct 4 wires. Probe's stem Ø3 mm, length 230 mm. 4-wire connection cable with connector, length 2 metres.
- **TP47.1000.0** Immersion probe, sensor Pt1000. Probe's stem Ø3 mm, length 230 mm. 4-wire connection cable with connector, length 2 metres.
- **TP87.100.0** Immersion probe, sensor Pt100 direct 4 wires. Probe's stem Ø3 mm, length 70 mm. 4-wire connection cable with connector, length 1 metres.
- **TP87.1000.0** Immersion probe, sensor Pt1000. Probe's stem Ø3 mm, length 70 mm. 4-wire connection cable with connector, length 1 metres.
- **TP47** Only connector for probe connection: Pt100 direct 4 wires, Pt1000 2 wires.

DELTA OHM metrology laboratories LAT N° 124 are ISO/IEC 17025 accredited by ACCREDIA for Temperature, Humidity, Pressure, Photometry / Radiometry, Acoustics and Air Velocity. They can supply calibration certificates for the accredited quantities.

# Notes

# WARRANTY

The manufacturer is required to respond to the "factory warranty" only in those cases provided by Legislative Decree 6 September 2005 - n. 206. Each instrument is sold after rigorous inspections; if any manufacturing defect is found, it is necessary to contact the distributor where the instrument was purchased from. During the warranty period (24 months from the date of invoice) any manufacturing defects found will be repaired free of charge. Misuse, wear, neglect, lack or inefficient maintenance as well as theft and damage during transport are excluded. Warranty does not apply if changes, tampering or unauthorized repairs are made on the product. Solutions, probes, electrodes and microphones are not guaranteed as the improper use, even for a few minutes, may cause irreparable damages.

The manufacturer repairs the products that show defects of construction in accordance with the terms and conditions of warranty included in the manual of the product. For any dispute, the competent court is the Court of Padua. The Italian law and the "Convention on Contracts for the International Sales of Goods" apply.

# **TECHNICAL INFORMATION**

The quality level of our instruments is the result of the continuous product development. This may lead to differences between the information reported in the manual and the instrument you have purchased.

We reserves the right to change technical specifications and dimensions to fit the product requirements without prior notice.

## **DISPOSAL INFORMATION**



Electrical and electronic equipment marked with specific symbol in compliance with 2012/19/EU Directive must be disposed of separately from household waste. European users can hand them over to the dealer or to the manufacturer when purchasing a new electrical and electronic equipment, or to a WEEE collection point designated by local authorities. Illegal disposal is punished by law.

Disposing of electrical and electronic equipment separately from normal waste helps to preserve natural resources and allows materials to be recycled in an environmentally friendly way without risks to human health.



CE RoHS